



X-ray Topography system based on Lang method

XRT-100

Operating Manual

Manual No. ME14903B01

Rigaku Corporation

Contents

2. Concept of X-ray diffraction topography and possible experimental setup	3
2.1 Lang method	4
2.2 Step scanning section topography.....	4
2.3 Surface reflection topography.....	5
3. Specifications.....	6
3.1 X-ray generator	6
3.2 Four-inch Lang camera	6
3.3 Controller and software	7
3.4 X-ray Vision Camera	8
3.5 Mount and radiation enclosure.....	8
4. Installation conditions	9
4.1 Ambient environment	9
4.2 Power supply	9
5. Structure and Functions	10
5.1 Incident optics.....	11
5.2 Topographic goniometer (topogonio).....	12
5.3 Sample holder	14
5.4 Positioning the film cassette bracket and film cassette.....	16
5.5 Receiving optics.....	17
5.6 X-ray Vision Camera	18
5.7 Controller and motor-driven axes	19
6. Operation	20
6.1 Starting up and shutting down the system	20
6.2 Lang method	21
6.2.1 Principle of Lang method.....	21
6.2.2 Principles of forming lattice defect images.....	22
6.2.3 Spatial resolution.....	25
6.2.4 Procedures.....	26
6.2.5 Simultaneous reflection image.....	31
6.2.6 Scattered X-ray shielding plate for transmission topography	33
6.3 Section topography	34
6.3.1 Principles of section topography	34
6.3.2 Tools for section topography	35
6.3.3 Procedures for section topography	35
6.4 Surface reflection topography.....	37
6.4.1 Principles of surface reflection topography	37
6.4.2 Procedures for surface reflection topography	39
6.4.3 Scattered X-ray shielding plate for reflection topography	42
6.5 Darkroom equipment	43

6.5.1	X-ray film and nuclear plates	43
6.5.2	Developer	43
6.5.3	Darkroom equipment	43
7.	Maintenance.....	45
7.1	Replacing the mylar films on the sample holder rings	45
7.2	Installation	46
Appendix	47
A	Data	47
B	Guide for references related to topography	49

1. Outline

Technological improvements in the electronics industry have made it possible to produce a variety of large single crystals, and such crystals are now being used in semiconductors and other devices. This unit is an X-ray diffraction topography system that enables you to nondestructively observe lattice defects such as dislocations, impurity segregation/precipitation, and stacking fault in such large crystals.

The basic features of this system are as follows:

- (1) It allows large samples (up to 4" wafers) to be studied;
- (2) There is no contamination of the sample by virtue of the Mylar-foil sample sandwich method;
- (3) There is no fear of X-ray exposure, as crystal alignment can be done with a remote ω -angle and in-plane rotation mechanism;
- (4) Alignment is easy with the X-ray Vision Camera and the vertically movable mechanism of the X-ray detector; and
- (5) Normal focus is available as an X-ray source.



Fig.1(a) Overview of XRT-100 system



Fig.1(b) XRT-100 main body

2. Concept of X-ray diffraction topography and possible experimental setup

X-ray topography is a measurement technique whereby lattice defects and strain in a single crystal are observed as an X-ray diffraction image. The acquired X-ray diffraction image is called an X-ray topograph. As is widely known, the sample crystal diffracts the incident X-ray beam when the crystal is positioned so that lattice planes with lattice spacing d satisfy the Bragg condition, $2d\sin\theta_B = \lambda$, with respect to wavelength λ of the incident X-rays. If the X-ray radiation field ABCD includes structural defects as shown in Fig. 2.1, the diffracted X-ray intensities differ between beams from areas containing defects and from areas free of defects. The difference is recorded on X-ray film or other media to produce an X-ray image corresponding in point to point fashion with the crystal. This image is the X-ray topograph.

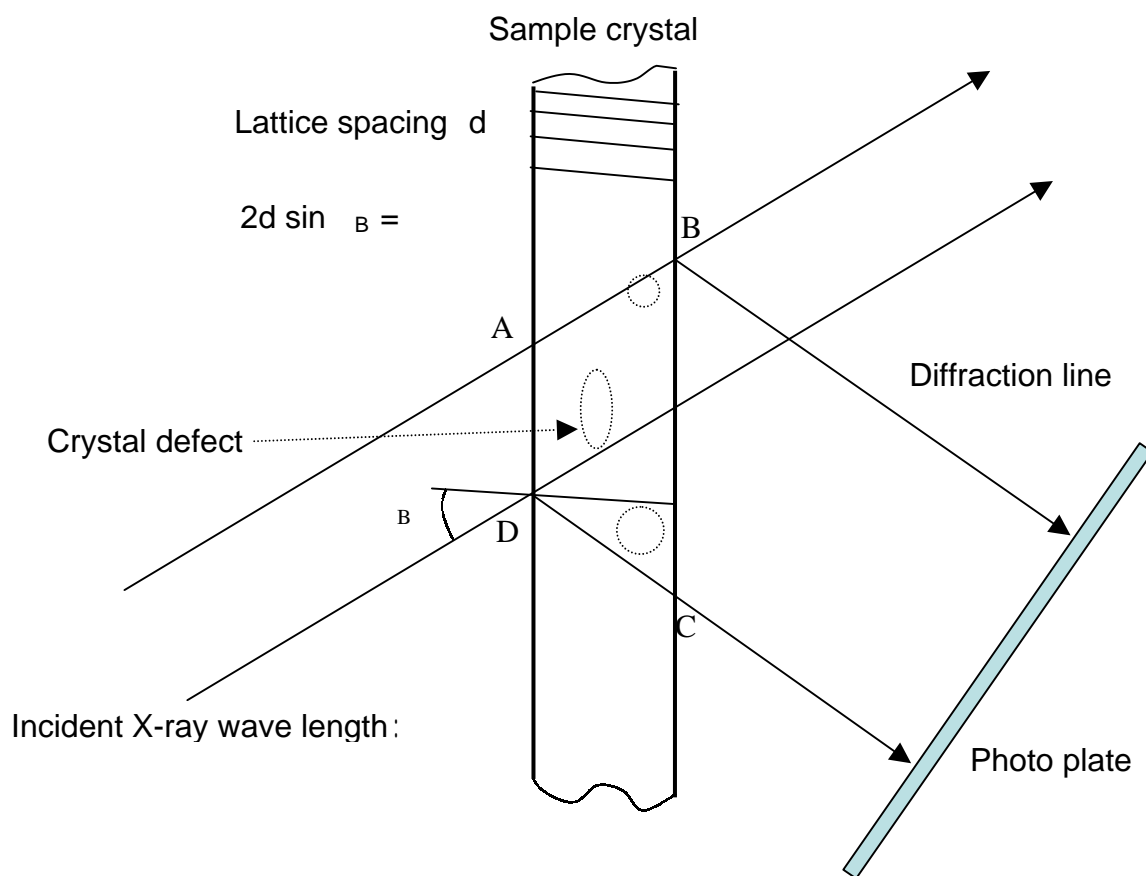


Fig.2.1 Concept of X-ray topography

The following subsections describe the topography techniques supported by the XRT-100 system.

2.1 Lang method

Lang method uses a Mo target and performs topography in transmission geometry. This method is capable of capturing an image of defect distributions throughout the wafer and can be applied for wafer quality inspections. Topography with a fine focus produces high spatial resolution.

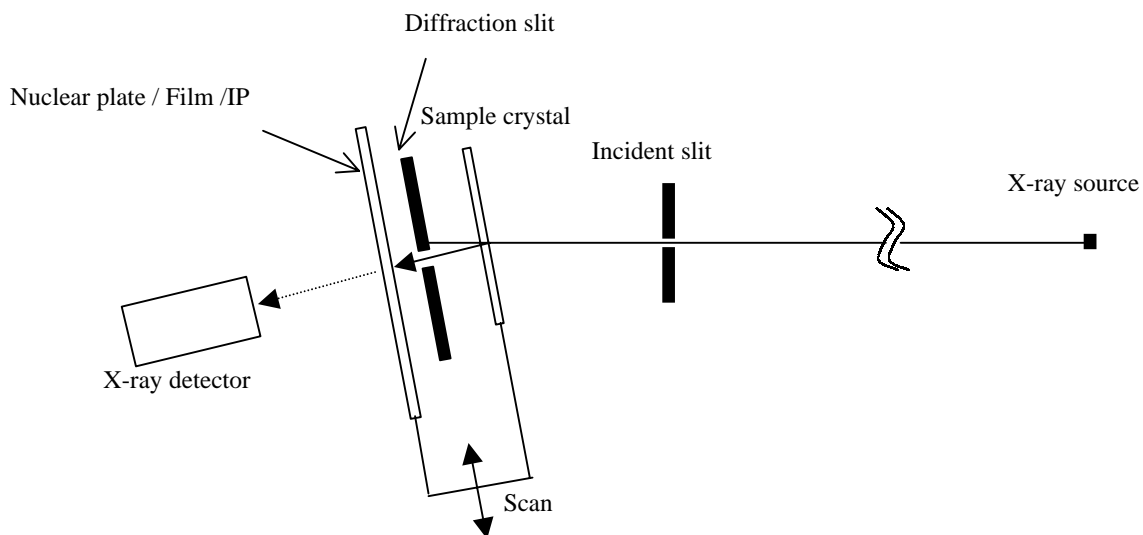


Fig. 2.2 Lang method

2.2 Step scanning section topography

Step scanning section topography uses a Mo target and performs topography in transmission geometry with a 10- μm slit placed near the sample, the sample, and the nuclear plate fixed for time exposures. This method provides information on defect distributions in the axis of thickness of the sample. The stage can be moved stepwise to make multiple exposures.

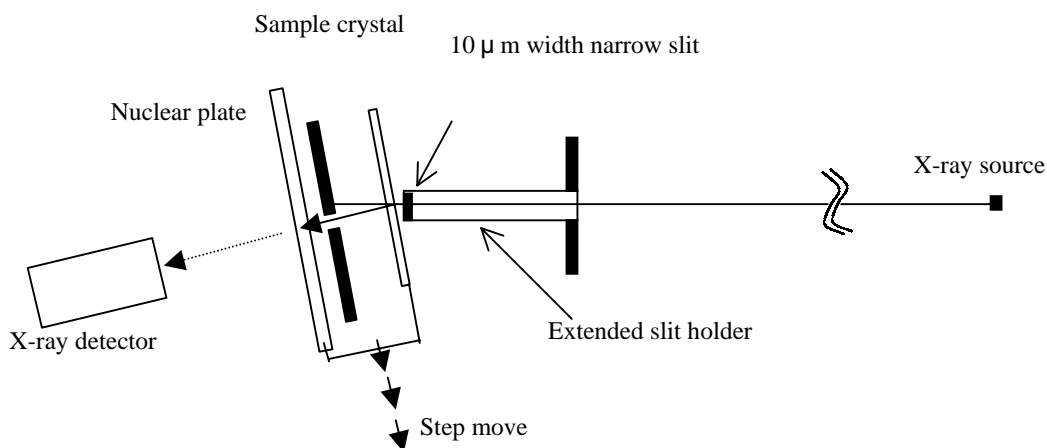


Fig. 2.3 Step scanning section topography

2.3 Surface reflection topography

Surface reflection topography uses a Cu target and performs topography in reflection geometry. This method is capable of imaging the entire wafer and can be used for crystal quality inspections. It provides information on surface defects.

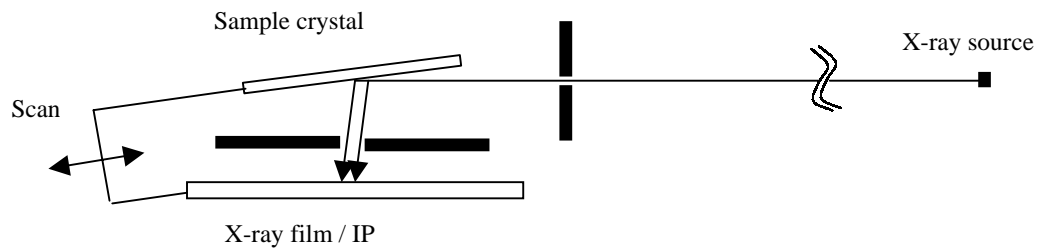


Fig. 2.4 Surface reflection topography

3. Specifications

3.1 X-ray generator

- (1) Maximum rated output: 3 kW
- (2) Rated tube voltage: 20 kV to 60 kV (1-kV step)
- (3) Rated tube current: 2 mA to 50 mA (1-mA step)
- (4) Stability (tube voltage and tube current): $\pm 0.01\%$ or less (power supply fluctuations should be $\pm 10\%$ or less.)
- (5) Tube shield: Horizontal
- (6) X-ray window: Equipped with electromagnetic shutter
- (7) Target (selected): Mo(for transmission topography) and Cu(for reflection topography) selected
- (8) Safety device: Cooling water error (water quantity, water pressure, water temperature)
XG overload fault (overload, line current, filament current)
Tube voltage HIGH/LOW fault
Filament burnout detection
Emergency stop switch
Ground-fault circuit breaker
- (9) Water circulation pump: Water circulation pump for 3-kW equipment.
Temperature control: $\pm 0.5^\circ\text{C}$

3.2 Four-inch Lang camera

- (1) Distance between X-ray source and sample: 1,100 mm
- (2) Center height of X-ray beam above mount: 490 mm
- (3) 2θ rotation plate Set range: 0° to 115° , can be set manually. Read with degree scale
- (4) ω rotation plate
Coarse adjustment: 0° to 150° can be set manually. Read with degree scale
Fine adjustment: Pulse motor operable within the range of $\pm 3.5^\circ$.
Resolution: 0.0001 deg/step
- (5) Sample holder: Usable both transmission and reflection topography
 ϕ rotation : In-plane rotation of 360° for both transmission and reflection topography. Pulse motor-driven.
Holder: The sample is inserted between mylar sheets.
Tilting: $\pm 5^\circ$, manual setting.
- (6) Scanning stage: Scanning range of ± 52 mm. Pulse motor-driven.

- (7) X-ray detector holder: Fixed on the 2θ rotation plate. The vertical positions of the X-ray detectors can be set by manual in a range of ± 50 mm from the height of the X-ray beam center.
- (8) First slit: 3 mm width. Fixed slit in the cylinder pipe.
- (9) Second slit
- (i) Width-limiting slit : Straight slits with widths of 0.1 mm, 0.2 mm, 0.5 mm, 1 mm, 1.5 mm, and 2 mm
- (ii) Height-limiting slit: 25 mm, 50 mm, and 75 mm
- (iii) Slit for section topography: Equipped with an extension slit holder to reduce the distance between the slit and the sample. Slit widths of 0.01 mm (for topography), 0.2 mm and 1 mm (for setting). Slit height of 20 mm.
- (iv) Slit for zero setting: 0.2 mm
- (10) Third slit: Width from 0.5 mm to 16 mm. Pulse motor-driven.
- (11) Film cassette: 4.5-inch square film cassette. 1-inch \times 1.5-inch nuclear plate cassette. Two sets are supplied.
- (12) Connection fitting: Connection fitting between the Lang camera and the X-ray tube shutter.

3.3 Controller and software

- (1) Control motor axes: Five axes, including ω fine adjustment, ϕ , scanning stage, third slit (right and left).
- (2) Curve correction: The scanning stage and the ω fine adjustment are moved together for curve corrections. The minimum curve correction radius is 5 m.
- (3) Communication method: RS232C
- (4) Topography control software:
- Axis setup
 - Traverse topography measurement
 - Curve correction data measurement
 - Section topography measurement
- (5) Personal computer section:
- HP desktop PC, d530
 - Pentium 4 2.8 GHz
 - 512 MB Memory
 - 80 GB HDD
 - 48 \times CD-RW
 - OS Windows 2000
 - 17-inch TFT Monitor
 - Color printer

3.4 X-ray Vision Camera

The X-ray Vision Camera mounted on the counter holder is used for sample axial adjustments and optical adjustments. X-ray images are transferred to the PC through a video capture board and viewed on the PC CRT monitor.

- (1) Configuration
 - X-ray image intensifier
 - CCD video camera
 - Video capture board
- (2) Sensitive area: Approximate diameter of 40 mm
- (3) Luminance gain: 12,000
- (4) Resolution limit: 100 μm
- (5) CCD: 2/3-inch

3.5 Mount and radiation enclosure

- | | |
|-------------------------------------|--|
| (1) Mount dimensions: | 1,000 mm W \times 1,000 mm D \times 720 mm H |
| (2) Radiation enclosure dimensions: | 1,400 mm W \times 1,425 mm D \times 2,000 mm |
| (3) Radiation enclosure system: | Iron plate radiation enclosure. Equipped with doors at the front and rear. Equipped with a lead plate measuring 3-mm in thickness (along the axis of the direct beam). |
| (4) Leakage X-ray: | 2.0 μSv or less |
| (5) Warning lamp: | Mounted on the top of the enclosure |

4. Installation conditions

4.1 Ambient environment

- (1) Temperature: Within the range from 15°C to 25°C.
- (2) Temperature fluctuation: Within the range of $\pm 0.5^\circ\text{C}$ or less during measurement. To ensure thermal stability of mechanical components, including the goniometer, temperature fluctuations should be kept as low as possible, even when no measurement is taking place.
- (3) Vibration: No perceptible vibrations.

4.2 Power supply

- (1) X-ray generator: Three-phase 200 VAC, 20 A
- (2) Controller and personal computer section: Single-phase 100 VAC, 15 A
- (3) X-ray Vision Camera: Branched from the power supply (3)
- (4) Air-cooled water circulation pump: Three-phase 200 VAC, 15 A
- (5) Grounding: Two series of Class-D grounding (One is used solely for the computer)

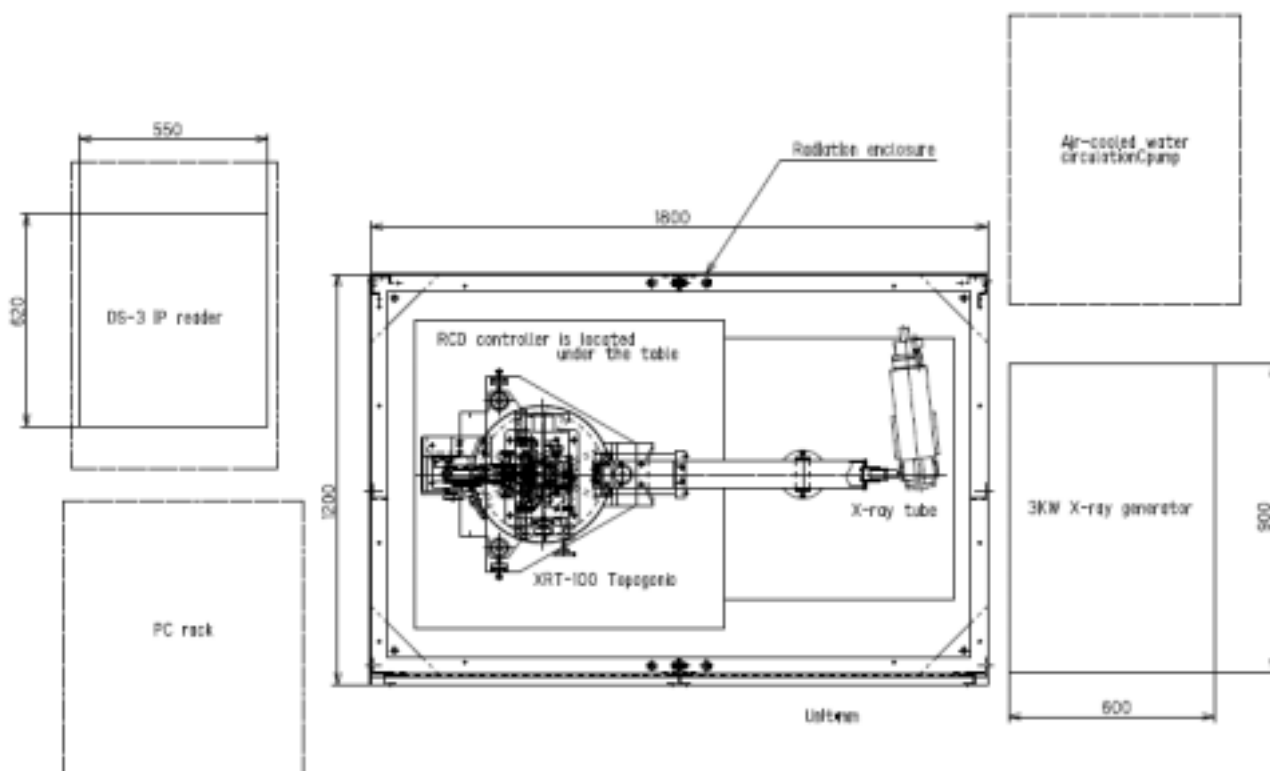


Fig. 4.1 Configuration of XRT-100 system

5. Structure and Functions

Fig. 5.1 provides a view of the entire XRT-100CCM system.

The equipment consists of the following five sections:

- (1) 3kW X-ray generator
- (2) Main unit
- (3) Controller and Control PC
- (4) X-ray vision camera

The X-ray generator and the main unit are each covered by a radiation enclosure. The main unit consists of the incident optics, topographic goniometer (topogoniometer), and the receiving optics.

The X-ray source is placed on the X-ray generator work table. The X-ray beam emitted from the X-ray source passes through the slit system to form a parallel beam and is directed to the sample attached to the topographic goniometer (topogoniometer). The diffracted X-rays are detected by the receiving optics, including the X-ray vision camera. The system can be controlled from the control PC.



Fig. 5.1 View of entire XRT-100CCM system

5.1 Incident optics

(1) X-ray pass and first slit

Set the Lang camera at a glancing angle of about 6° relative to the target face, then couple the X-ray pass with the X-ray window through the protector (Fig.5.2). The X-ray pass is a shield cylinder. Turn the shield cylinder so that the clamp dial of the protector directs upward. The first slit is fixed in the shield cylinder. The slit width is 3 mm. The purpose of the first slit is to keep excessive X-ray radiation from reaching the camera.

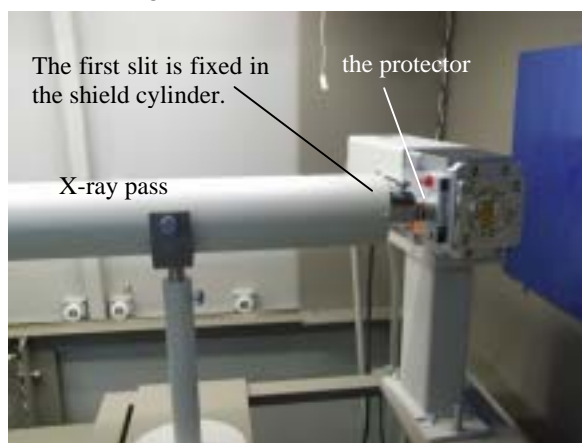


Fig. 5.2 X-ray pass and connecting section

(2) Second slit, S2

The second slit is used to restrict the radiation field of the X-ray beam incident on the crystal. Use a width-limiting slit and a height-limiting slit based on the size of the crystal. Insert the slits into the slit box.

Besides these slits, slits for section topography and a setting slit are provided.



Fig. 5.3 Second slit section

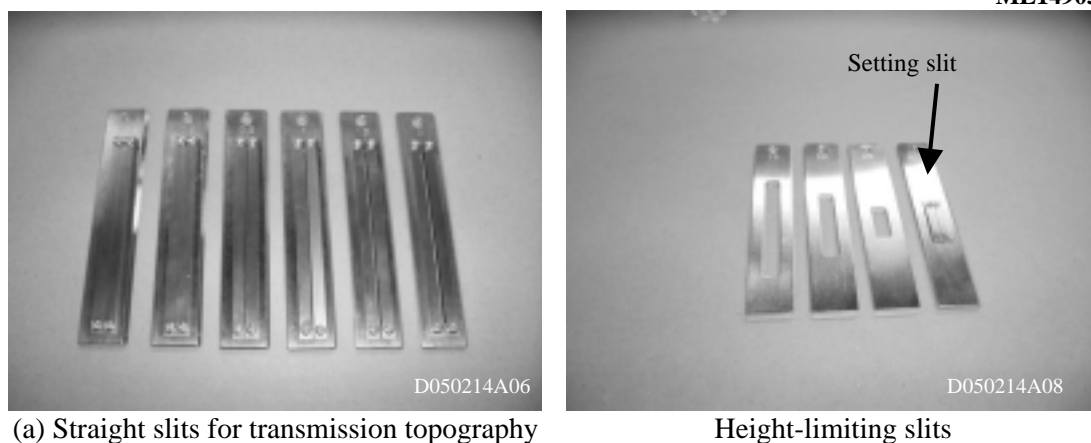


Fig. 5.4 2nd Slits

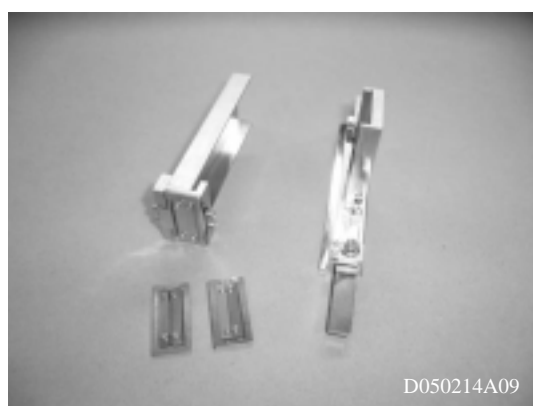


Fig. 5.5 Slits for section topography

5.2 Topographic goniometer (topogonio)

(1) 2θ rotation plate

The 2θ rotation plate is a rotation mechanism that holds the X-ray detectors and the third slit. It rotates around the topogonio center and sets the detectors and the slit to the diffraction angle (2θ). Set the angle with the scale plate and the reference line.

(2) ω rotation plate

The ω rotation plate holds the scanning stage and the sample holder and rotates around the topogonio center. The plate is a rotation mechanism for adjusting the incident angle of the X-ray beam with respect to the sample crystal. Loosen the clamp screw to rotate the plate manually. Make coarse settings for the desired angular position with the scale plate and the reference line. Tighten the clamp screw to enable fine adjustment by the tangent bar method. A single increment of the ω vernier scale corresponds to an angle of 0.01° .

(3) Scanning stage

The scanning stage is a precision sliding stage for acquiring wide-area topographs. It holds the crystal and the film and slides back and forth along an axis parallel to the crystal plane. Its range of motion is ± 52 mm. If you try to move the stage beyond this limit, an over-run microswitch (limiter) will activate and halt movement.

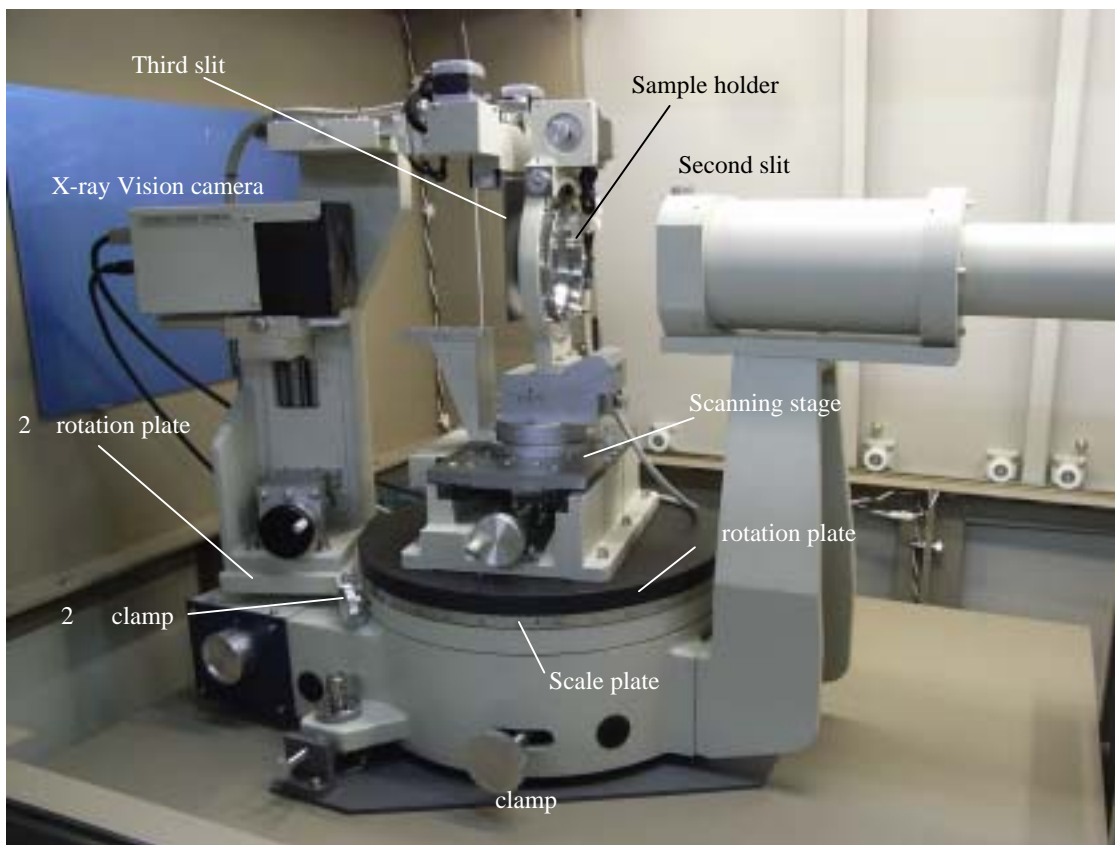


Fig. 5.6 Topographic goniometer (topogoniometer)

5.3 Sample holder

Fig. 5.9 is a photograph of the sample holder section. The sample holder is intended to hold a wafer or slab sample measuring less than 1 mm thick. The sample holder employed the Mylar-foil sample sandwich method.

The sample can be held between the Mylar-foil that is extended on a pair of holder rings (Fig .5.7). The holder rings can be used for both transmission and reflection topography.



Fig. 5.7 Sample holder ring

Set the crystal into the holder rings as described below.

- (a) Place the crystal wafer in the center of Ring A (Fig. 5.8 (a)). Crystal thickness must be less than 1 mm.
- (b) Place Ring B (Fig. 5.8 (b)) over Ring A and press. The crystal is held in place by the spring effect.
- (c) Insert Rings A and B (Fig. 5.8 (c)) into the sample holder. Insert them until they stop; at this position, the crystal plane will match the rotation axis of the goniometer. The spring effect holds the crystal at an arbitrary in-plane rotational position.
- (d) Fig. 5.81 (c) indicates the relationship between the X-ray incidence from the front and rear of the crystal.

The sample holder has a mechanism for adjusting the crystal orientation, permitting both in-plane rotation and tilting.

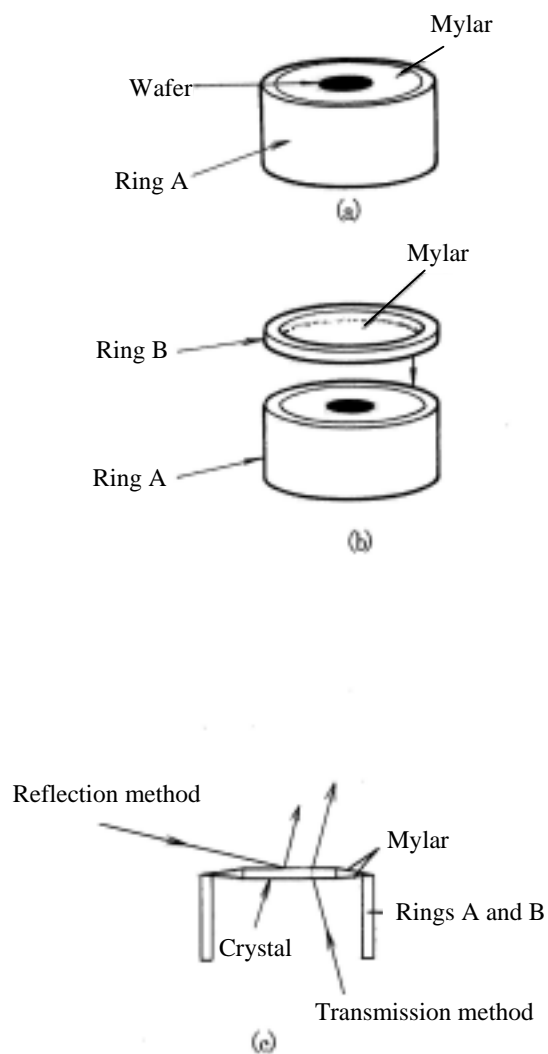


Fig. 5.8 Sample holder rings



Fig. 5.9 Sample holder with holder rings inserted

5.4 Positioning the film cassette bracket and film cassette

The film cassette is held by the cassette bracket via the film cassette holder. The cassette bracket can be rotated around the sample. This rotation is called ψ .

Set the film cassette parallel to the sample and the surface of the third slit to shorten the distance between the sample and the film.

Set the cassette bracket angle to zero. The sample holder features a scale. Set the ψ angle to $\psi = 0$. The film cassette holder becomes parallel to the sample and the surface of the third slit. Move the cassette a little. This moving records the topograph in the center of the film. The amount of movement is calculated by the equation $Lc \cdot \tan(2 -)$. Lc is a camera length that is a distance between the sample and the film.

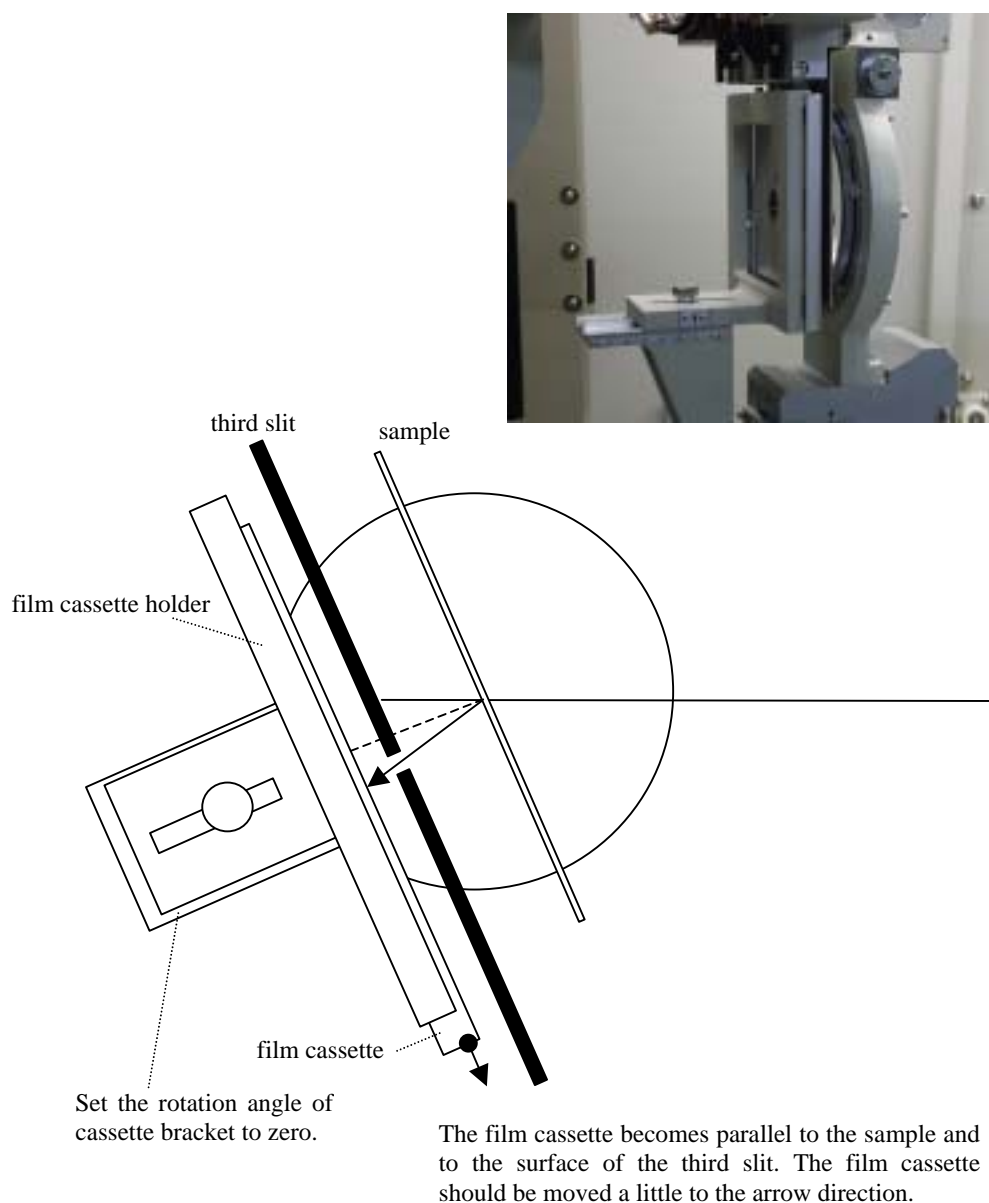


Fig. 5.10 Positioning the film cassette bracket and film cassette

5.5 Receiving optics

(1) Third slit (variable slit) (Fig. 5.11)

The third slit is used to reduce the effects of beam scattering on the film and to provide protection against direct exposure to X-ray beams. The width of the slit is independently variable, driven by a pulse motor. The entire slit can be set rotated from the position perpendicular to the diffracted beam within the range of approximately $\pm 30^\circ$. This mechanism makes it possible to set the scanning direction of the scanning stage parallel to the surface of the third slit and to reduce the distance between the sample and the film cassette. The film cassette and width of the third slit will be large when the Lang camera is used with a large sample; this means the resulting distance between the sample and the film is large, frequently resulting in poor resolution. The rotational setting of the third slit may help remedy this shortcoming.



Fig. 5.11 Third slit

(2) X-ray detector and movement mechanism

We used the X-ray vision camera as an X-ray detector. You can move the sensitive area of the X-ray vision camera and set vertically to maximize the sensitive area. This is a convenient way to make axial adjustments of the large crystal in topography geometry.

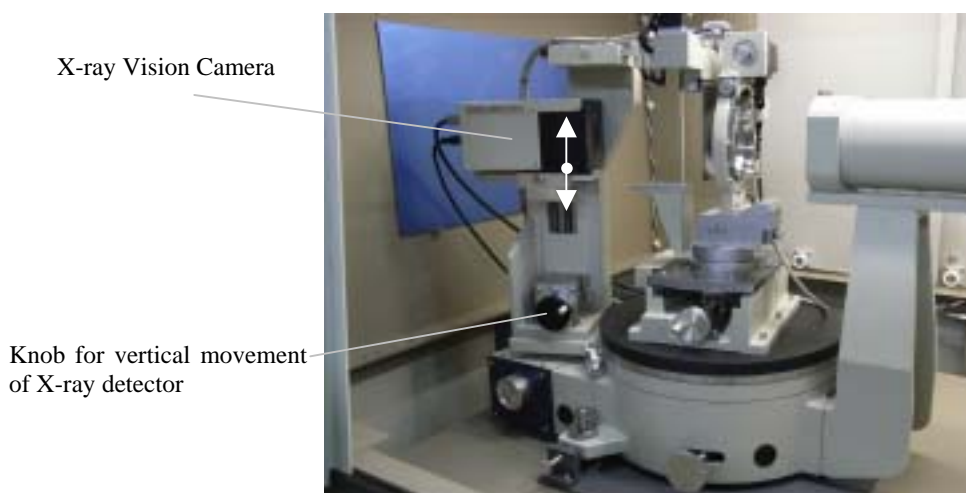


Fig. 5.12 X-ray detector and movement mechanism

5.6 X-ray Vision Camera

The X-ray Vision Camera is a highly sensitive X-ray TV that converts a faint X-ray diffraction image into visible light on a fluorescent screen, amplifies the image with the image intensifier (I.I.), and captures the output image with the lens-coupled CCD. The CCD operates at the TV rate. The output signal can be observed on the video monitor. Operate the topographic goniometer remotely while observing the video monitor and use the camera to search for diffraction lines and to perform axial adjustments of the crystal and other adjustments.

Operate the X-ray Vision Camera as follows: Turn on controller power. Set the Gain control to zero. High voltage is supplied to the I.I. only after this step. Turn the Gain control clockwise to increase sensitivity. For standard use, set the Gain to 5 on the scale. Adjust the value according to X-ray intensity.

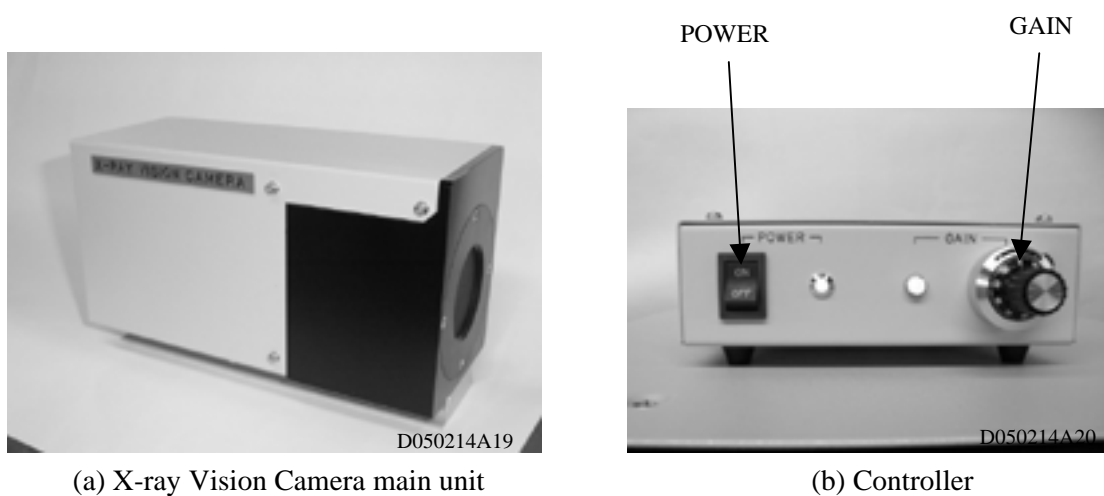


Fig. 5.13 X-ray Vision Camera

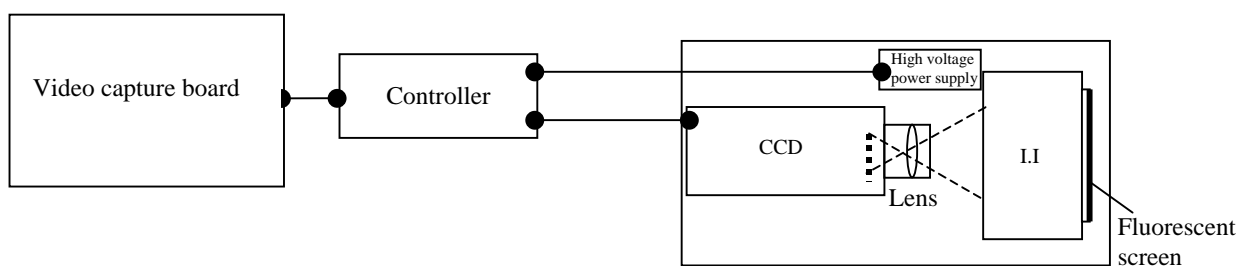


Fig. 5.14 Block diagram of X-ray Vision Camera

5.7 Controller and motor-driven axes

The XRT-100 system uses the RCD3, Rigaku's standard controller. The following five axes are motor-driven: ω , stage, ϕ , variable slit left and variable slit right. Fig. 5.20 schematically illustrates the arrangement of the motor-driven axes with respect to the incident X-ray direction when the 2θ and ω axes are set to zero. The arrows indicate the positive direction of translation or rotation.

Table 5.1 shows the range of movement and resolution of the axes.

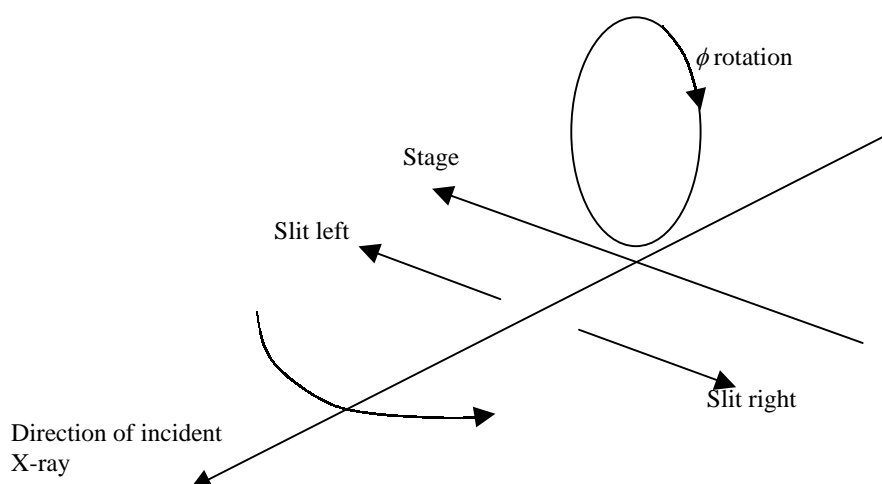


Fig. 5.15 Configuration of motor-driven axes and positive direction of translation and rotation (indicated by arrows)

Table 5.1 List of motor-driven axes

Name of axis	Range of movement	Resolution
ω axis	$\pm 3.5^\circ$	0.0001°
Stage	± 52 mm	0.001 mm
ϕ axis	$\pm 360^\circ$	0.005°
Variable slit left	-5 ~ +8 mm	0.00005 mm
Variable slit right	-5 ~ +8 mm	0.00005 mm

6. Operation

This chapter describes the procedures for operating the XRT-100 system, starting with procedures related to daily startup and shutdown. The various topographic imaging techniques are described thereafter.

For discussions of general equipment handling, refer to the Lang Camera for Large Samples XRT-100 software Instruction Manual (ME14902E01).

6.1 Starting up and shutting down the system

Start up the system as follows:

- (1) Provide power to the system. Turn on the breakers for the water circulation pump (three-phase, 200 V), X-ray generator (three-phase, 200 V), and the personal computer and controller (single-phase, 100 V).
- (2) Turn on power for the TV camera system. Set the Gain of the X-ray Vision Camera controller to zero, then to around 5 on the scale.
- (3) Start the PC and PC peripheral devices. (Power on in the following order: printer, monitor, PC.) Launch the topography control software, and begin initialization.
- (5) Power on the X-ray generator. The water circulation pump will begin operating, and the system will move to the Ready state. (The doors of the radiation enclosure must be closed for Ready state.)
- (7) Turn on X-ray generation. Warm up the equipment for approximately 5 minutes at minimum load.
- (8) Gradually increase the X-ray load voltage (kV) and current (mA) manually. Take time to set the load to the value desired for measurement.
- (9) For operations, launch the topography control software.
Shut down the system as follows (generally speaking, shutdown procedures are simply the startup procedures reversed):
 - (1) Decrease the X-ray load, starting with current (mA), to the minimum load of 20 kV and 2 mA.
 - (2) Turn off X-ray generation. Allow the system to stand for approximately 3 minutes until the cooling water has automatically stopped.
 - (3) Turn off power for the X-ray generator.
 - (4) Close the topography control software and turn off power for the PC and peripheral devices.
 - (5) Turn off the power switch for the TV camera system.
 - (6) Turn off the 100-V power supply, then turn off power for the X-ray generator and the water circulation pump.

6.2 Lang method

6.2.1 Principle of Lang method

Lang method, the topography method currently in widest use, was developed by A.R. Lang. Fig. 6.1 illustrates the underlying principles. The X-rays emitted from a point X-ray source are directed and width-limited by a long second slit and allowed to strike the sample, irradiating the sample crystal along a narrow strip. When the in-plane orientation and the incident angle are adjusted to satisfy the diffraction conditions for the lattice plane of the crystal, diffraction occurs across the entire area irradiated.

An X-ray tube with an Ag, Mo, or Cu cathode is generally used as the X-ray source.

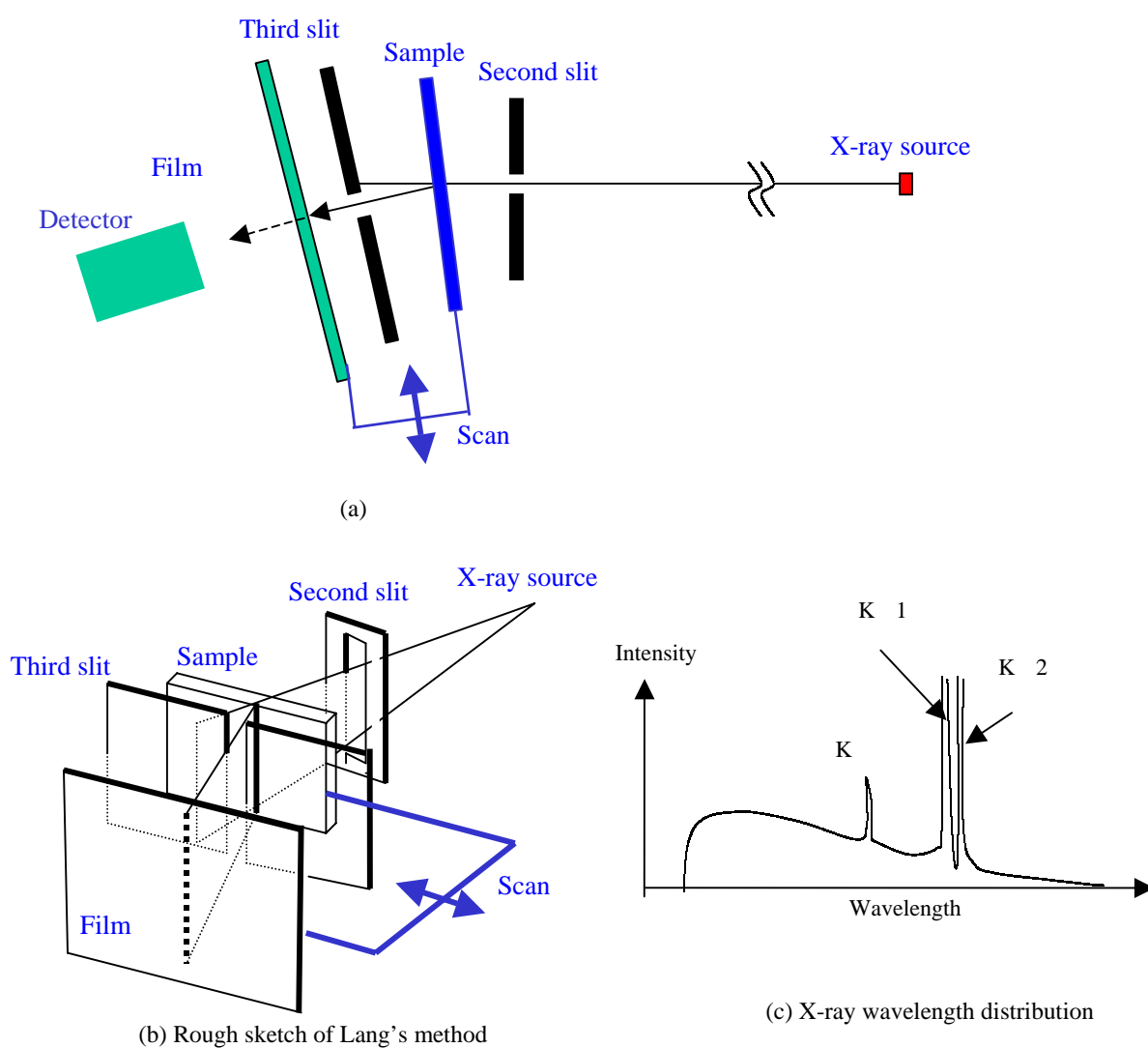


Fig. 6.1 Principle of Lang's method

The third slit has a function that reduces the background generated by the scattered X-rays. Narrow the width of the slit so that it shields the direct X-ray beam through the sample and allows only the diffracted X-ray beam to pass.

Scan the sample crystal parallel to the sample surface together with the film in the above configuration to obtain a diffraction image over the entire sample.

When capturing a topograph, place the film as close to the sample as possible to produce a high-resolution image. A topograph acquired in this manner is known as a traverse topograph. Since it projects a three-dimensional defect image to a two-dimensional image, it is also called a projection topograph.

6.2.2 Principles of forming lattice defect images

(1) Extinction contrast

When the incident X-ray beam is perfectly monochromatic and collimated, the angular width of the diffracted X-ray beam from a perfect crystal will be several seconds or less. The incident X-ray beam used with Lang's method has an adequately large divergence angle of several minutes. Among incident X-rays, those satisfying the precise diffraction conditions excite the transmission and diffraction waves and are reflected from the lattice planes repeatedly, as shown in Fig. 6.2 (a).

The standing wave formed by two waves with the same amplitude propagates parallel to the lattice plane from the incident point O to B. As the waves shift from the diffraction conditions, the transmitted and diffracted waves acquire different amplitudes and spread in the triangle OAC (referred to as the Borrmann fan) formed by the axis of transmission OA and axis of diffraction OC.

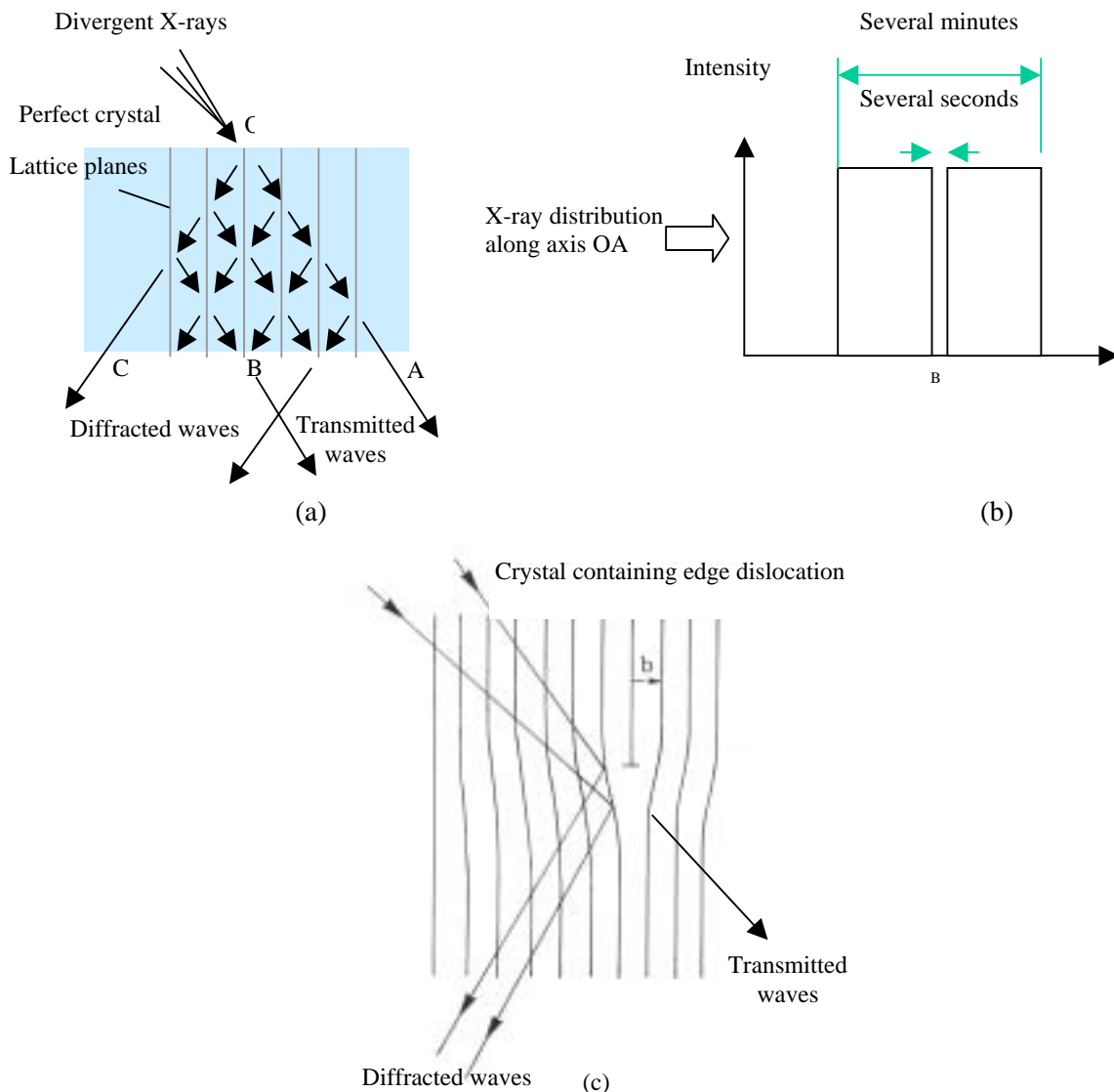


Fig. 6.2 Extinction contrast

Only a portion of the incident X-rays propagate between A and C; most X-rays propagate along the axis of transmission OA, as shown in Fig. 6.2 (b). The integrated diffraction intensity from a perfect crystal is weak.

However, as shown in Fig. 6.2(c), if the crystal contains a dislocation, the lattice planes are bent by several seconds with a change in the lattice spacing in the region within several $10\ \mu\text{m}$ of the dislocation core. When the crystal is translated and the dislocation crosses the X-rays along the axis OA, the X-rays incident at various axes satisfy the diffraction conditions around the dislocation and are diffracted. The diffracted X-rays diverge, and the integrated intensity becomes extremely large. This is why the Lang topograph contains a strong image of the dislocation over the weak image of the perfect portion of the crystal with weak diffraction intensity.

The phenomenon whereby diffraction intensity is weaker for a crystal with few flaws is known as the extinction effect. The image formation principle is called the extinction contrast. The extinction contrast is formed by a thin crystal satisfying the condition $\mu t < 1$, where μ is the linear absorption coefficient and t the thickness.



Fig. 6.3 Example of extinction contrast (positive) topography of dislocations in Si crystal

(2) Anomalous transmission contrast

For thick samples for which $\mu t \gg 1$, anomalous transmission contrast is formed. Here, if the crystal is perfect, the X-rays that propagate along axis OB, in Fig. 6.2 (a) experience only an anomalously small absorption and pass through and reach the rear side of the crystal. X-rays propagating in other directions are almost entirely absorbed. If the crystal contains a dislocation along axis OB, the multiple scattering stops, and X-rays do not pass through the crystal. Thus, the topograph exhibits a contrast pattern that is the reverse of the extinction contrast pattern. Fig. 6.4 illustrates this behavior schematically.

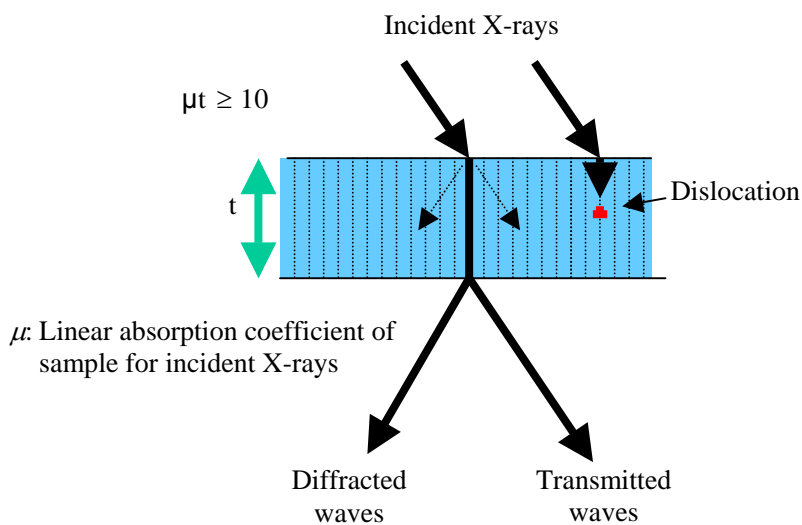


Fig. 6.4 Anomalous transmission contrast

Fig. 6.5 is an X-ray topograph of a GaAs (100) 2-inch wafer doped with In, showing images of the dislocation and slip bands formed by anomalous transmission contrast.

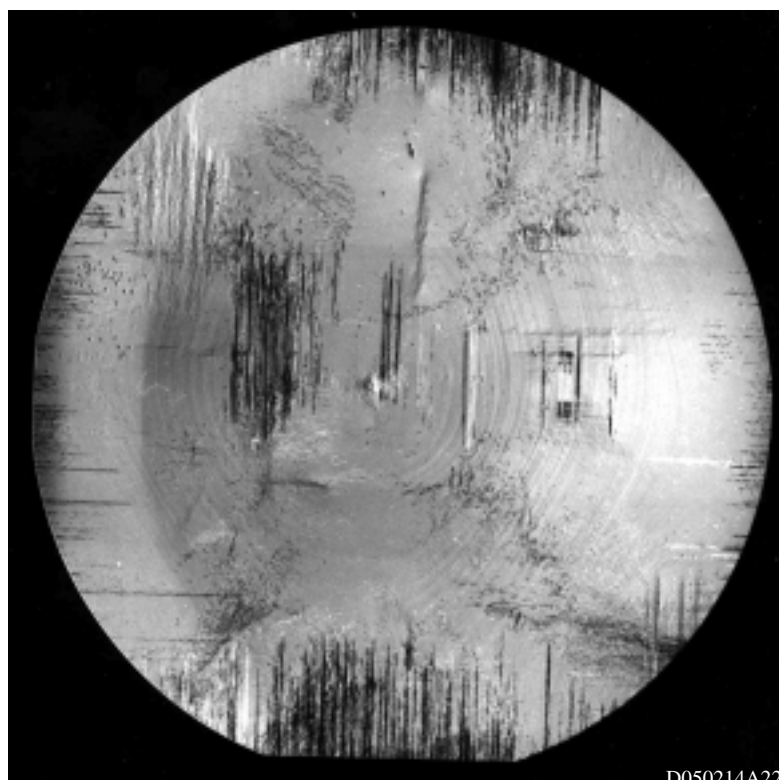


Fig. 6.5 Example of anomalous transmission contrast (positive)

Orientation contrast is another contrast type. If the crystal has several domains differing in crystal orientation or contains a lineage (an imperfect structure wherein the orientation of the crystal changes continuously along a linear boundary), the degree of the diffraction condition satisfied changes, changing the resulting intensity.

6.2.3 Spatial resolution

The magnification ratio for X-ray topography is basically 1. Resolution depends on the size of the silver halide grains of the X-ray films used. Nuclear emulsion plates are used to record high-resolution topographs. For the Type L4 plates made by Ilford, the size of the silver halide grains is approximately $0.2 \mu\text{m}$ after development. Topographs made using these plates can be magnified by a factor of around 100.

Fig. 6.6 illustrates geometric spatial resolution. For the sake of simplicity, the diffraction angle (2θ) is assumed to be zero. The resolution in the dimension of height, R_h , is a simple geometric unsharpness and is given by $R_h = F_h \cdot m/l$. Here, F_h is the height of the X-ray source, l the distance between the X-ray source and the sample, and m the distance between the sample and the film. The resolution in the dimension of width, R_w , is given by $R_w = m \cdot \Delta\theta$, where the divergence angle from the defect is $\Delta\theta$. The divergence angle is dominated by angular spreading due to wavelength dispersion and is given by $\Delta\theta = (\Delta\lambda/\lambda)\tan\theta_B$. Here, $\Delta\lambda$ is the natural width of the characteristic X-rays and θ_B the Bragg angle. $l \cdot \Delta\theta$ is generally smaller than width F_w of the X-ray source, and the resolution in the dimension of width is independent of the size of the X-ray source. To optimize the system setting for high resolution, use an X-ray source having a micro focus of $50 \mu\text{m}$ and take the parameters at approximately $l = 500 \text{ mm}$ and $m = 10 \text{ mm}$. This achieves a resolution of $1 \mu\text{m}$ to $2 \mu\text{m}$ along both height and width.

The XRT-100 system is designed for large aperture measurements and uses a strong normal-focus X-ray generator having a normal focus of 0.5 mm for F_h , to balance against exposure time. Here, $l = 1,100 \text{ mm}$ or greater and approximately $m = 20 \text{ mm}$, and the resolution in the dimension of height degrades to approximately $10 \mu\text{m}$. In practice, this should not present a problem.

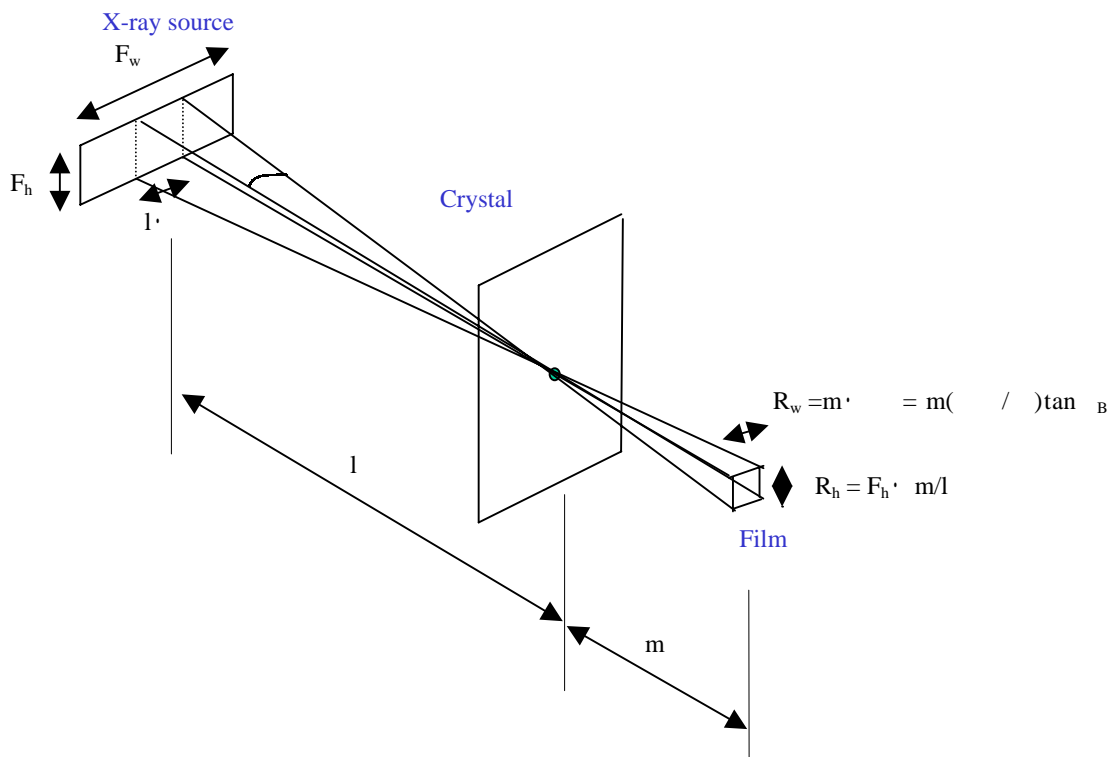


Fig. 6.6 Geometric resolution

6.2.4 Procedures

This section describes the procedures for various steps ranging from attaching the sample to measuring topography. For information on using the personal computer, refer to the Lang camera for Large Samples XRT-100 software Instruction Manual (ME14902E01).

(1) Attaching the sample

Determine the orientation of the sample in advance (e.g., by the Laue method). When measuring an industrial silicon wafer, the orientation flat (OF) indicates the in-plane orientation.

Hold the sample in the sample holder rings by following the producers described in Section 5.3. Adjust the in-plane orientation so that the intended diffraction plane is nearly vertical. Press the holder ring into the sample holder until it stops. The sample surface will be parallel to the scanning direction and incorporate the axis of the ω rotation plate.

(2) Preparations

Set the 2θ rotation plate to the predetermined 2θ angle. Set the vertical position of the X-ray Vision Camera to the zero position.

The width of the second slit should be approximately equal to the apparent width of the X-ray focus (normally 1 mm). Use a height-limiting slit suitable for the size of the sample. Open the third slit wide. Remove the film cassette and film cassette holder and set the scanning table to the center position.

Make coarse settings to set the ω rotation plate at roughly the angular position satisfying the diffraction conditions. (Clamp the stage firmly with the clamping screw).

For transmission topography, as the diffraction plane, select a lattice plane nearly perpendicular to the sample surface. Set the ω angle as $\omega = \theta_B$ (Bragg angle) if the diffraction plane is perpendicular to the sample surface. Denoting the angle of tilt from the sample surface normal as δ , set the ω angle as $\omega = \theta_B \pm \delta$. The sign depends on δ , the direction of the angle of tilt.

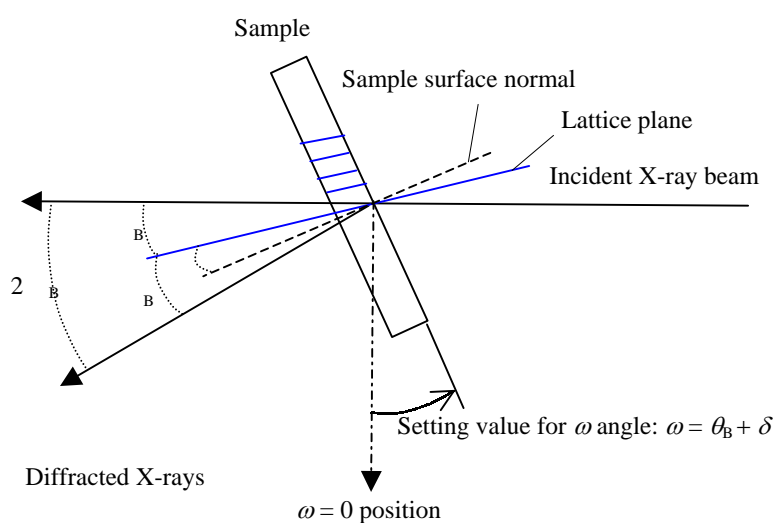


Fig. 6.7 Setting value for ω angle

As an example, Table 6.1 lists the calculated results for the possible transmission geometry for a Si wafer of (001) plane orientation. The values are for MoK α_1 ($\lambda = 0.070926$ nm) X-rays.

Table 6.1 Reflection indices suitable for transmission topography of Si wafer with (001) plane orientation

hkl	d - (nm)	Unit: Degree			
		2			
111	0.313556	6.49	12.99	35.24	41.76
220	0.19201	10.64	21.29	0	10.65
311	0.16375	12.51	25.02	17.55	30.06
400	0.13577	15.14	30.28	0	15.14
331	0.12459	16.54	33.07	13.26	29.80
422	0.11086	18.66	37.31	24.10	42.75
511	0.10452	19.83	39.67	11.10	30.93
333	0.10452	19.83	39.67	35.26	55.10
531	0.09180	22.72	45.45	9.73	32.46
602	0.08586	24.39	48.78	18.43	42.82


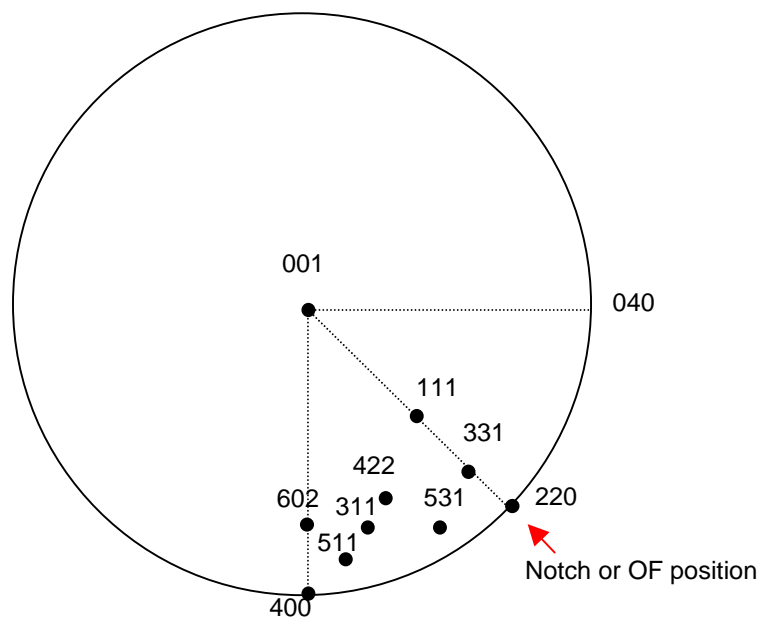


Fig. 6.8 Stereogram for (001) wafer


(3) Axial adjustment for sample crystal

Generate X-rays and perform axial adjustments for the crystal. To do this, start “Axis set up” (PC operation) and operate the system while observing the X-ray diffraction image with the X-ray Vision Camera. Suppress X-ray output so that the output signal from the X-ray Vision Camera does not saturate. The diffracted X-ray intensity will depend on the sample and the diffraction vector selected. Appropriate X-ray output adjustments will require some operator experience.

Open the X-ray shutter, perform the ω scan, and search for a diffracted beam. Fig. 6.9 (a) shows the output image just after the image of the diffracted X-ray beams is captured. Since axial adjustments are incomplete, the diffracted beams in the image for the $K\alpha_1$ and $K\alpha_2$ appear as tilted lines. Set the orientation of the crystal by adjusting the in-plane rotation and by finely adjusting the ω angle so that the image is uniform from top to bottom, as shown in Fig. 6.9 (b). The field of view of the X-ray Vision Camera in the dimension of height is approximately 36 mm. If you are measuring a large sample, move the detector vertically while observing the image of the diffracted beam to confirm that the diffraction condition is achieved across the full vertical range of the sample. Make fine orientation adjustments, if necessary.

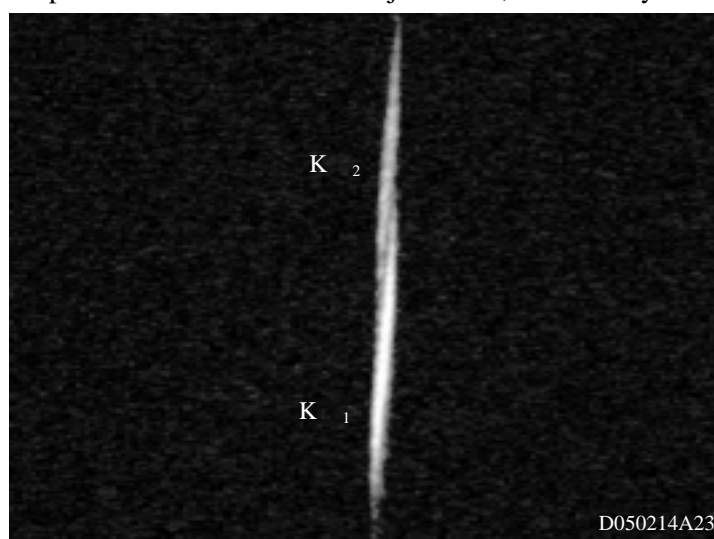

(a) Image of diffracted beam captured by ω scan

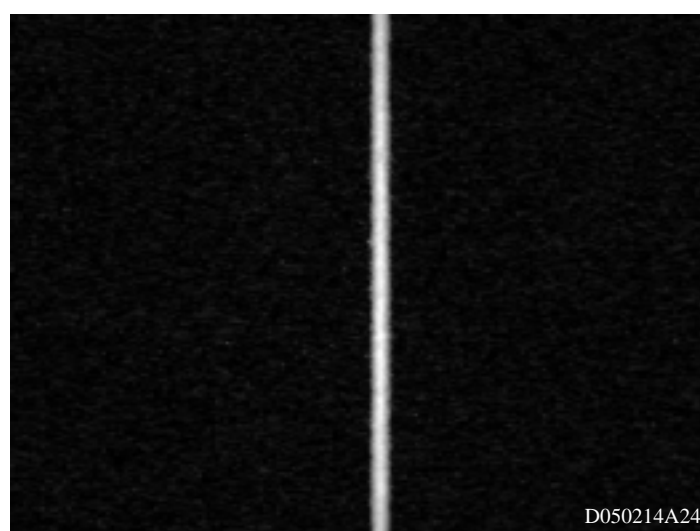

(b) Image after adjusting orientation

Fig. 6.9 Axial adjustment

- (4) Determining the position of the third slit
Close the X-ray shutter and move the third slit closer to the crystal. Adjust the setting so that the crystal surface and the slit surface are more or less parallel. Position them as close as possible while avoiding collision between the crystal and the slit as the crystal moves back and forth.
- (5) Setting the width of the third slit
Open the X-ray shutter once again and observe the image of the diffracted X-ray beam. Select the adjustment axis from “slit right” and “slit left,” moving the slits to narrow the slit. After confirming that the slits block the image of the diffracted X-ray beam, set the slits to make the slit width slightly larger than at the position where the image is blocked. Adjust the position of the other slits in a similar manner.
- (6) Setting the scanning range
Move the scanning stage while observing the image of the diffracted X-ray beam. Continue moving the stage until the image of the diffracted X-ray beam disappears at the crystal edge. Stop the stage here and register the position of the crystal edge. If the stage is in the negative range, use the left edge register button. If the stage is in the positive range, use the right edge register button. When the crystal edge positions are registered, the XRT-100 system will automatically update scanning range settings for traverse topography. If the diffracted X-ray beam shifts from the diffraction condition as the scanning table moves, the sample is warped. In this case, follow the instructions for acquiring the curve correction data given under Item (7).
- (7) Setting sample curve correction data
If the sample is bent, set the sample curve correction data. By following the procedures indicated in the operation program move the sample position in stepwise increments, acquire ω value for the precise Bragg condition and crystal edge information by ω scan adjustments, and register the information. Use the acquired step data to interpolate the data. Select the interpolation method from the spline function and first-degree to fifth-degree polynomial interpolation based on the least mean squares method. When the first-degree polynomial is selected, the average curvature radius is calculated and displayed.
If the sample is warped significantly, ω correction may fail in the exposure step for the traverse topography described in Item (9), and the equipment may generate an error and stop. Curve corrections are possible for samples with a curvature radius of 5 m or greater.
- (8) Positioning the film cassette (Fig. 6.10)
Insert an X-ray film or nuclear emulsion plate into the film cassette and mount on the camera. Insert the film cassette into the cassette holder, support with the cassette bracket, and fasten in place with the clamp screw. Set the film cassette as close to the sample as possible. For this purpose, set the film cassette as close to perfectly parallel to the surface of the third slit as possible while avoiding collisions with the third slit within the scanning range for the scanning stage. Read the distance between the sample and the film with the reference lines and scale.



Fig. 6.10 Positioning the film cassette

(9) Exposure

Set the X-ray output to the maximum value. Set the exposure conditions and begin the exposure. The exposure time will depend on exposure conditions: scanning range, scanning speed, and number of scans. Since the diffraction intensity depends on the X-ray load, sample, and diffraction vector and since exposure time depends on the sensitivity of the X-ray film or nuclear emulsion plate, establishing suitable exposure conditions often requires some user experience.

(10) Development

Use the specified developer and fixer solutions to develop the X-ray film or nuclear emulsion plate. Develop X-ray films in the developer at 20°C for 5 minutes, rinse in water or soak in a 1.5% glacial acetic acid solution (stop bath) for approximately 30 seconds, then fix in a fixer solution at 20°C for 5 minutes. Finally, rinse the film in water for approximately 30 minutes.

Developing a single side of the X-ray film will facilitate magnified observation. X-ray films typically have structures in which a photographic emulsion slightly thicker than 10 μm is applied to both sides of a polyester film base. If the image is magnified after development of both sides, the resulting image will appear in double. To prevent this, develop only one side of the film. Use a wide adhesive tape (such as Scotch Book Tape 845 or vinyl tape) to cover the rear side of the film, then proceed with development. After soaking the film in the stop bath, remove the tape and fix the film.

Proceed with care when developing a nuclear emulsion plate with a thick emulsion layer. If you attempt to develop the plate in a warm environment, the emulsion may peel off. To prevent this, maintain the temperature of the developer solution at 4°C, place the plate in the solution, and develop for approximately 1 hour. Following development, rinse the plate with water for approximately 30 seconds, then thoroughly fix the plate in fresh fixer solution at 20°C. Rinse the plate with water for approximately 1 hour. Dry the rinsed plate in air. Note that accelerated drying may create cracks in or cause the emulsion to peel from the glass substrate.

6.2.5 Simultaneous reflection image

A Lang topograph will sometimes contain streaks along the dimension of width. This is a diffraction phenomenon known as “simultaneous reflection” which always appears when the sample has a large aperture.

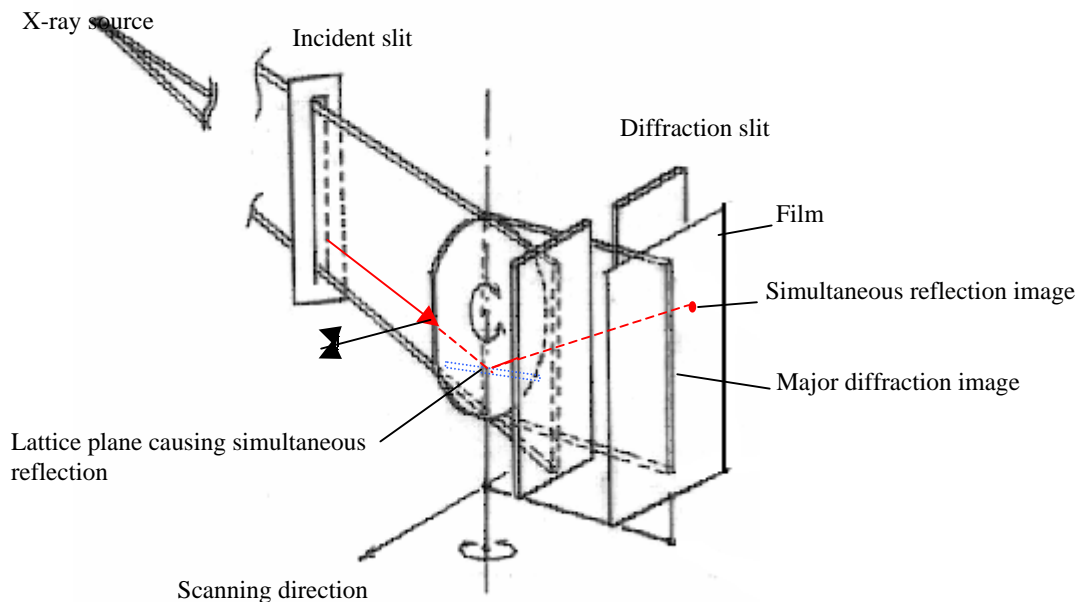


Fig. 6.11 Rough sketch of simultaneous reflection

As shown in Fig. 6.11, X-rays emitted from the X-ray source are shaped by the long incident slit and irradiated onto the sample. The ω angle and ϕ angle are adjusted to set the orientation of the sample so that the diffraction conditions on the specified lattice plane are satisfied. More specifically, the image of the diffracted X-ray beam is acquired under conditions whereby the lattice plane and longitudinal direction of the incident slit are parallel and the incident angle agrees with the Bragg angle. The image of the diffracted X-ray beam satisfies the diffraction condition across the full vertical range of the sample and is projected onto the film as a reed-shaped image.

In addition to this major diffraction image, the incident X-rays (arrow in the Fig.) may satisfy the diffraction condition for part of the sample. The diffracted beam is projected through the diffraction slit as a spot onto the X-ray film. This diffraction image occurs simultaneously with the primary image of the diffracted X-ray beam and is called a simultaneous reflection. A simultaneous reflection occurs because many lattice planes generating diffractions are oriented in several dimensions and directions in the crystal and because the incident beam diverges vertically. Simultaneous reflections are not restricted to characteristic X-rays; they can also occur with the wavelength component (Laue image) selected by the lattice plane causing the diffraction.

With Lang’s method, the sample and X-ray film are scanned parallel to the sample surface during exposure. The X-ray image corresponding to the outer shape of the sample is acquired by the primary image of the diffracted beam, and the simultaneous reflection image forms streaks along the dimension of width (scanning direction) that overlaps the primary image.

While the simultaneous reflection image overlaps the primary image of the diffracted beam, you will sometimes observe a white image. This is a residue of the simultaneous reflection.

Is it possible to prevent simultaneous reflections? Simultaneous reflections are basically a diffraction phenomenon and cannot be removed. However, one can move the position at which they appear via

rotation (χ rotation) around the normal of the lattice plane, as shown in Fig. 6.11. Setting strict (narrow) widths for the diffraction slit also helps reduce the number of the simultaneous reflections. The distance between the sample and film is called the camera length. Using a large camera length including the diffraction slit position can remove simultaneous reflections. Simultaneous reflections do not occur along the horizontal dimension; they are emitted downward or upward. Thus, the problem can be solved by increasing the camera length to reach a position at which a simultaneous reflection does not overlap the primary image of the diffracted beam. To optimize resolution, make the camera length as short as possible; long camera lengths are generally undesirable. It should be taken into consideration that simultaneous reflections are inevitable in high-resolution topography using X-ray film. However, since the resolution of topography with imaging plates (IP) is limited due to characteristics of IP, it may in certain cases be advantageous to increase the camera length to eliminate simultaneous reflections.

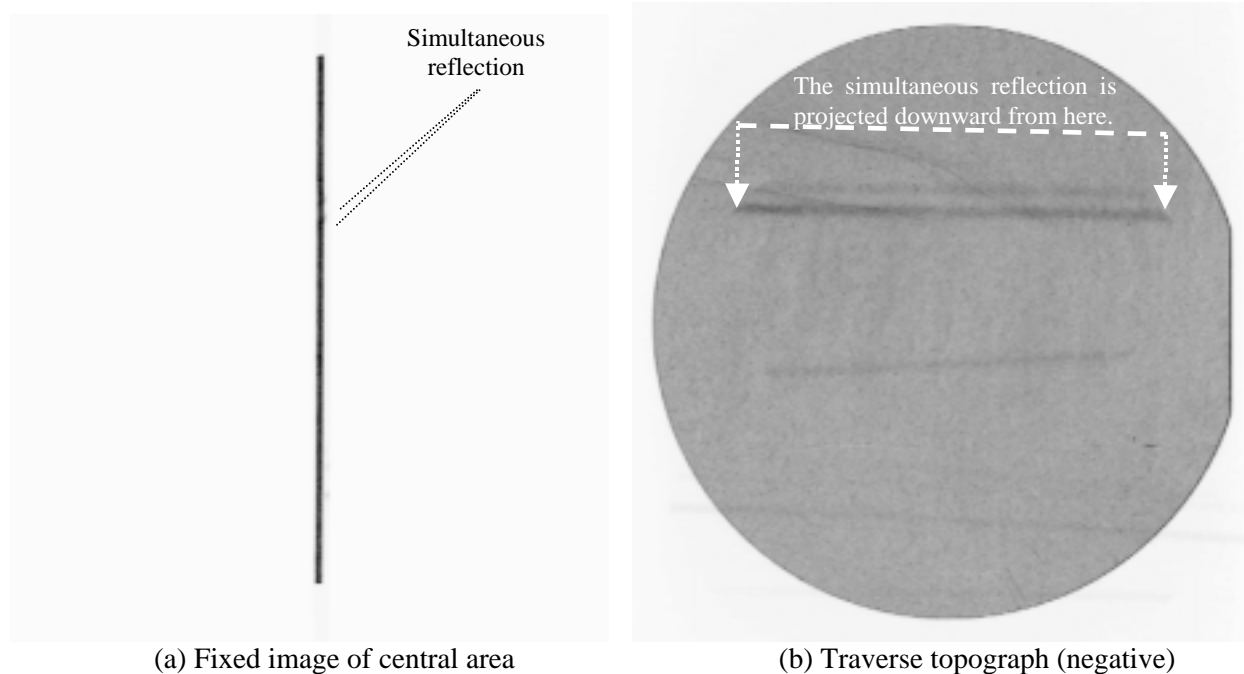


Fig. 6.12 Example of simultaneous reflection: (001) sapphire wafer with diffraction vector of 2-10

6.2.6 Scattered X-ray shielding plate for transmission topography

For transmission topography, attach the scattered X-ray shielding plate to obtain a clear topograph with low background. This method is useful when long exposure times are required for a thick sample or for crystals characterized by large absorption.

As shown in Fig. 6.13, incident X-rays are scattered by air or at the slit edges, generating scattered X-rays. These scattered X-rays pass through the third slit, reach the film, and are recorded as background. Use the scattered X-ray shielding plate to prevent this. The scattered X-ray shielding plate blocks scattered X-rays with its protruding wings and keeps them from reaching the film. Fig. 6.14 illustrates the effectiveness of using the scattered X-ray shielding plate, showing the background of a topograph captured without a scattered X-ray shielding plate.


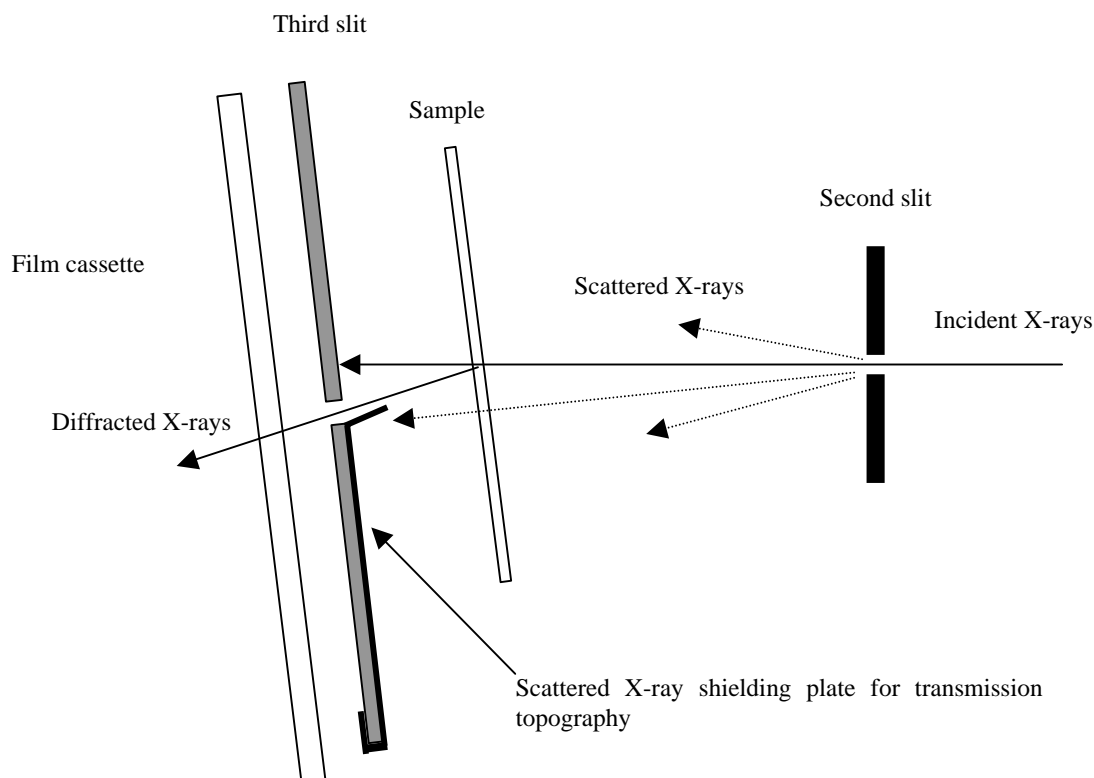


Fig. 6.13 Attaching the scattered X-ray shielding plate for transmission topography

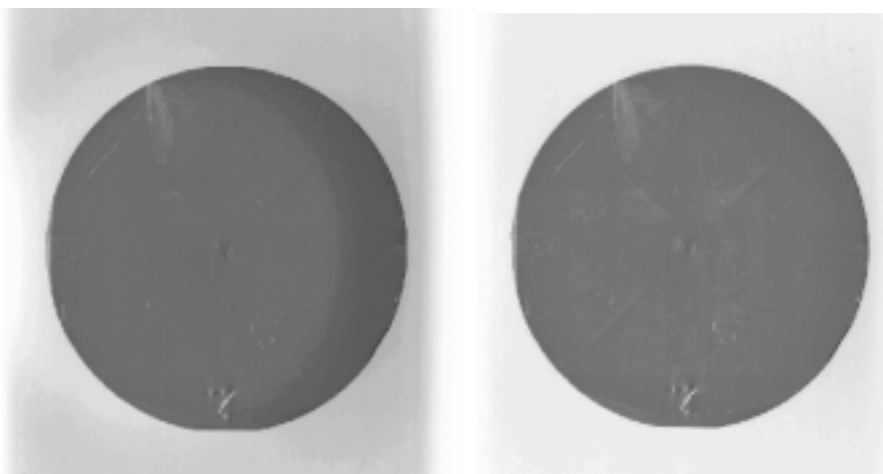

Without scattered X-ray shielding plate (D040318A04) With scattered X-ray shielding plate (D040318A05)

Fig. 6.14 Effectiveness of scattered X-ray shielding plate: GaAs $g = 220$ (negative)

6.3 Section topography

Section topography is a method for investigating the distribution of lattice defects along the dimension of depth by removing simultaneous reflections in a sample crystal. This section discusses section topography principles, tools, and imaging procedures.

6.3.1 Principles of section topography

X-ray diffraction image captures made with a stationary sample and stationary nuclear emulsion plate are referred to as section topography. In this case, the width of the incident X-rays is narrowed to approximately $10\ \mu\text{m}$ by the incident slit, which is placed close to the sample. A diffraction image captured with this method is called a section topograph.

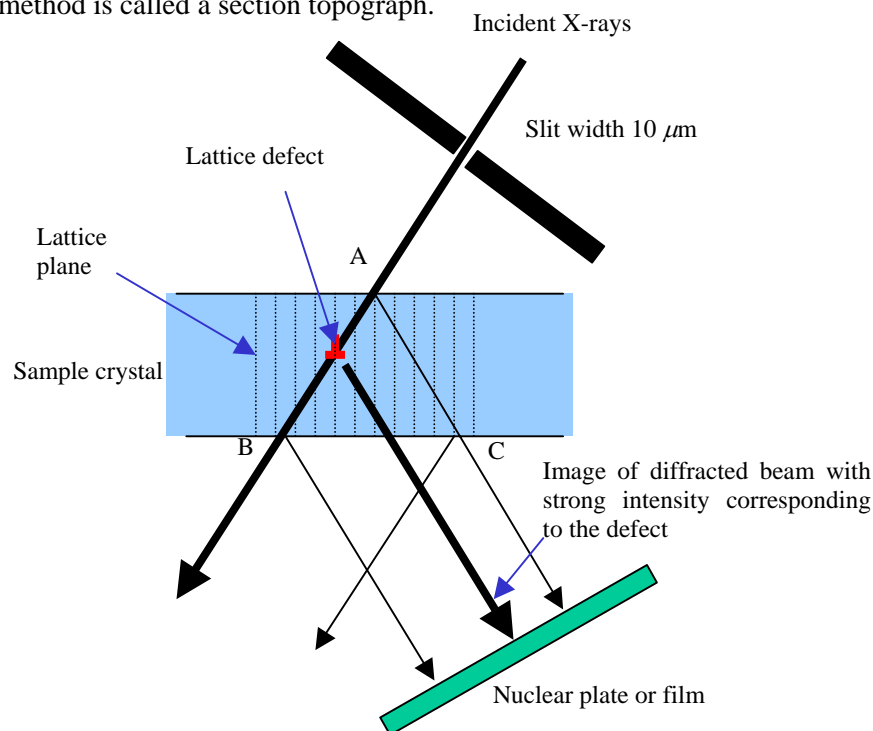


Fig. 6.15 Configuration of section topography (plan view)

Fig. 6.15 shows the configuration of a section topography in plan view. The X-ray incident beam is narrowed. If the diffraction conditions are met for the characteristic X-rays at the lattice plane of the sample crystal, the X-rays are diffracted. As the diffracted X-rays undergo multiple scattering in the crystal, they spread in the triangular region *ABC* and appear as an image with widths in the transmission and diffraction axes. If the crystal has relatively few flaws, interference fringes (Pendellösung fringes) will appear, caused by interference between the diffracted and transmitted waves. The visual appearance of the interference fringes is used to evaluate the perfection of the crystal.

When the crystal contains a dislocation in the direction of the direct X-ray beam as shown in the figure 6.15, a diffracted beam is generated nearby with strong intensity corresponding to the defect (defect image). The position of the defect image in the diffracted beam image recorded on the photographic plate provides information on the position of the defect with respect to depth. That is, it provides

information on the lattice defect distribution along the cross section of the crystal. Here, the term “section” is used to mean “cross section.”

In contrast to section topography, the ordinary topography measured by scanning is called traverse topography. Since the image is a projection of a three-dimensional image onto two dimensions, traverse topography is sometimes called projection topography.


6.3.2 Tools for section topography

To obtain a clear section topograph, note the following points:


- (i) Use a photographic plate with a fine-grain emulsion capable of producing high resolution (e.g., nuclear emulsion plate) (required).
- (ii) Use a narrow incident slit with a width of approximately 10 μm .
- (iii) Set the sample close to the 10- μm slit.

(If the distance between the sample and the slit is large, the angular divergence of the X-rays will diffuse the irradiated width even if the slit width is small, rendering the resulting section topograph unsharp.)

The XRT-100CCM system provides an extension slit holder for setting the sample close to the slit. The extension slit holder consists of a support attached to the second slit section and a slit holder into which the slit is inserted. The two components are secured by magnetism. The slits (setting slits) provided are 0.2 mm and 1 mm wide. Use these slits to create the Bragg condition and replace the slit into the 10- μm slit to capture the image. Select the 0.2 mm or the 1 mm slit as follows: If the X-ray source provides a normal focus, use the 1-mm slit; if it provides a fine focus, use the 0.2-mm slit.

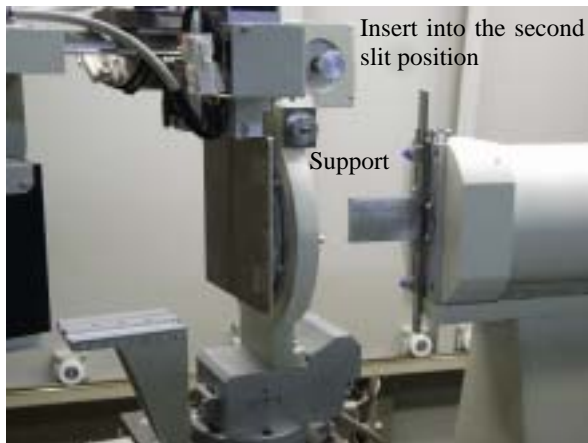

(a) Support

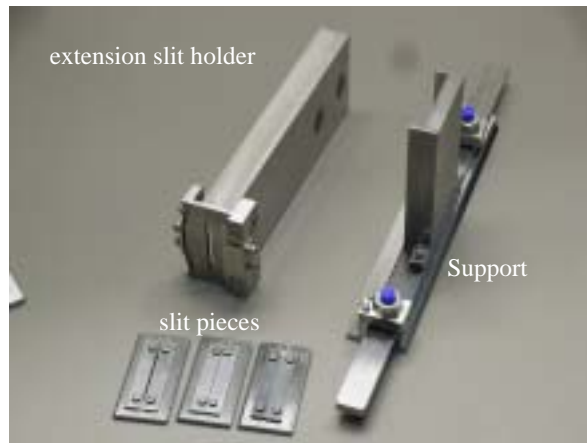

(b) Slit holder and slits

Fig. 6.16 Tools for section topography

6.3.3 Procedures for section topography

Before starting measurement, attach the sample, perform axial adjustments, and set the third slit as for ordinary transmission topography. Then, follow the subsequent procedures given below.

- (1) Attach the support for the extension slit to the second slit holder. Note that the support slides vertically. After attaching it, slide the support to set the appropriate vertical position.
- (2) Attach the setting slit to the slit holder. If the X-ray focus is normal, use the 1-mm slit. If it is fine

focus, use the 0.2-mm slit. Attach the slit to the support and move the slit close to the sample (Fig. 6.17).


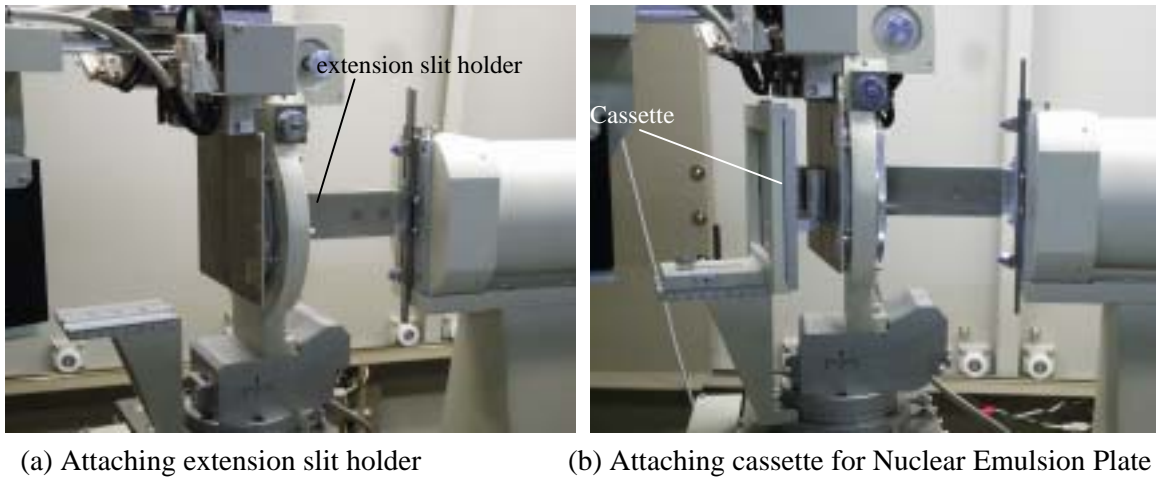


Fig. 6.17 Setting of section topography

- (3) Open the X-ray shutter and adjust the ω axis to the exact Bragg condition.
- (4) Replace the slits from the setting slit to the 0.01-mm (10- μ m) slit for section topography.
- (5) Attach the cassette and launch the control software for section topography and perform topography.

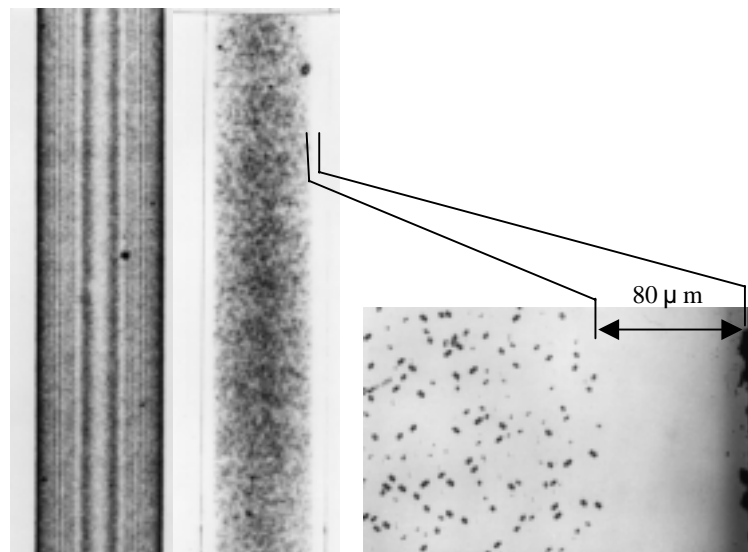

(a) Before heat treatment (b) After heat treatment (c) Optical microscopic image of etching pits in the cross section of the crystal after heat treatment
Section topograph

Fig. 6.18 Examples of section topograph (negative)


Fig. 6.18 (a) and (b) show section topographs of a Si-CZ crystal before and after heat treatment. Before heat treatment, we see X-ray interference fringes (Pendellösung fringes), indicating that the crystal is relatively free of flaws. Heat treatment precipitates SiO_x within the crystal, deforming the lattice and creating a granular contrast. As we see in the topograph, a denuded zone (D.Z.) is formed on the surface of the crystal by the out-diffusion of oxygen.

6.4 Surface reflection topography


6.4.1 Principles of surface reflection topography

Lang method (previously described in Section 6.2) is topography in the transmission geometry. Also possible is topography in the reflection geometry. Topography in the reflection geometry was first attempted by Berg in 1931 and refined by Barrett. Berg-Barrett topography is now virtually synonymous with reflection topography.

Fig. 6.19 shows a plan view and rough sketch of reflection topography. The X-rays emitted from a point X-ray source pass through the slit and irradiate the sample. Select a lattice plane that satisfies the diffraction condition with a small angle of incidence on the sample surface and a diffraction angle (2θ) close to 90 degrees. This setting makes adjustments easier. This type of reflection is known as asymmetric reflection. Asymmetric reflections spread the width of the image of the diffracted beams to approximately ten times the width of the X-ray source, enabling topography of a wide area of the crystal surface. Asymmetric reflections make it possible to set the sample close to the film, increasing resolution. It also makes it easier to position the receiving slit between the sample and the film. Scanning the sample in a manner similar to Lang method allows topography of the entire sample surface.

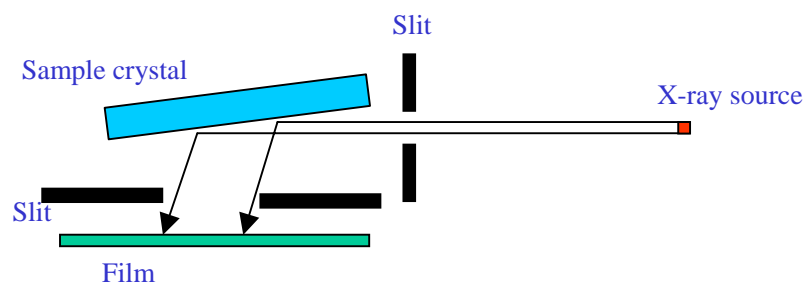

(a) Plan view

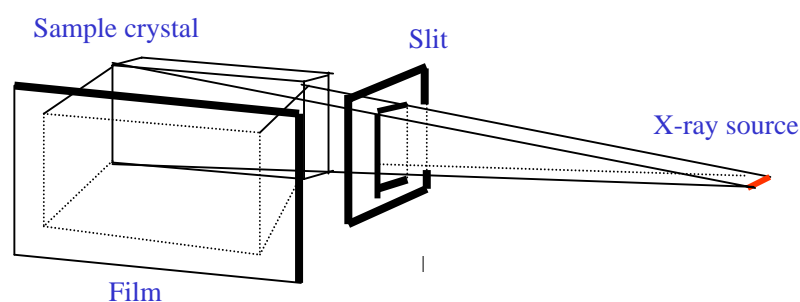

(b) Rough sketch

Fig. 6.19 Surface reflection topography

Berg-Barrett topography is free of any constraints on sample thickness and can be applied to examine a thick sample through which X-rays cannot penetrate. As with Lang's method, the principle of dislocation image formation is extinction contrast. In general, as the state of the crystal approaches perfection, diffraction intensity declines (a phenomenon known as the extinction effect). If such crystals contain lattice distortion due to lattice defects or other factors, the defective area generates intense diffracted X-ray beams, producing the contrast of the defect image and generating images of

dislocations on the surface of the crystal. However, the contrast of the dislocation image is weaker than with Lang method performed in transmission geometry.

If the crystal contains a region having a different orientation, this region has a different degree of diffraction condition satisfied and causes the intensity to change. This is a principle of image formation known as orientation contrast. It is suitable for observing a lineage, which is an assembly of dislocations, with orientation shifting in subgrains. Fluorite crystals often consist of subgrains with slightly different orientations. With such samples, orientation contrast makes it possible to observe the domain structure.

The method is also useful for the topography of heteroepitaxial films (thin film crystals deposited on a substrate crystal, with the same orientation but composition different from that of the substrate). The information acquired by surface reflection topography includes defects in layers within several to 100 μm from the surface.

Tables 6.2 to 6.4 show the reflection indices suitable for reflection topography for a Si wafer, a GaAs wafer and a fluorite crystal. The diffraction angle values are for the $\text{CuK}\alpha_1$ line.

The third slit and the film cassette have a rotation mechanism along the axis of the diffracted X-ray beam and a precession mechanism. Adjusting the rotation and precession reduces distortion in the topography due to the cosine effect.

Table 6.2 Reflection indices suitable for reflection topography of Si wafer with (001) plane orientation

Unit: degree

Reflection index	Bragg angle θ	Diffraction angle 2θ	Lattice plane tilting δ	Incident angle α
224	44.01	88.03	35.26	8.75
115	47.48	94.95	15.79	31.68
404	53.35	106.71	45	8.35

Table 6.3 Reflection indices suitable for reflection topography of GaAs wafer with (001) plane orientation

Unit: degree

Reflection index	Bragg angle θ	Diffraction angle 2θ	Lattice plane tilting δ	Incident angle α
224	41.87	83.74	35.26	6.6
115	45.07	90.13	15.8	29.27
404	50.42	100.83	45	5.41

Table 6.4 Reflection indices suitable for reflection topography of fluorite crystal with (111) plane orientation

Unit: degree

Reflection index	Bragg angle θ	Diffraction angle 2θ	Lattice plane tilting δ	Incident angle α
331	37.94	75.87	29.5	8.4
422	43.7	87.39	19.5	24.2
511	47.1	94.24	39.94	8.2
440	52.9	105.8	35.3	17.6

6.4.2 Procedures for surface reflection topography

This section describes procedures for reflection topography. Use the holder rings to grasp the sample, as in transmission topography.

(1) Considering reflection indices and setting the topography geometry

For samples whose crystal orientation is clear, prepare an angle table similar to Tables 6.1 to 6.3. This permits efficient setting of diffraction and incident and the in-plane rotational position. For samples whose crystal orientation is unknown, begin by measuring the orientation. Identify tilting angle δ of the lattice plane in question. If this is less than the Bragg angle, we can apply reflection diffraction geometry. The incident angle α is given by $\alpha = \theta_B - \delta$.

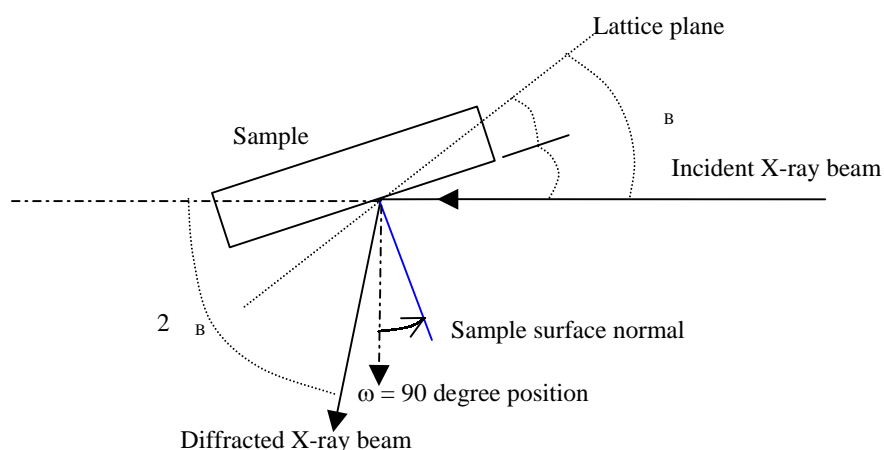



Fig. 6.20 ω angle setting for reflection geometry

Set the ω angle exceeding 90 degrees by further rotating it by the α angle: $\omega = 90 + \alpha = 90 + \theta_B - \delta$. Select the lattice plane for which the incident angle is less than 20 degrees to facilitate the setting of the topography geometry.


(2) Searching for reflection and axial adjustments

Set the diffraction angle 2θ to enable observations of the diffracted X-rays with the X-ray vision camera. To do this, set the lateral movement of the detector to the zero position.

Open the X-ray shutter, adjust the incident angle ω , and capture the diffracted beam. Adjust the in-plane rotation angle ϕ and the incident angle ω in alternation to extend the image of the diffracted beam vertically.

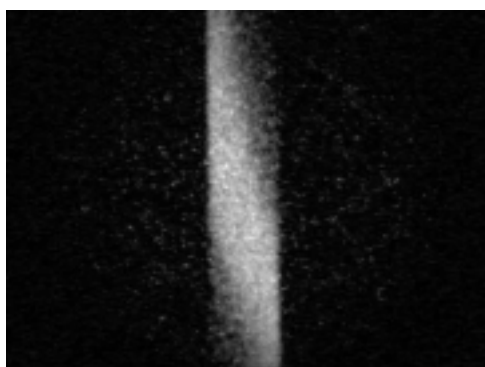

(a) Image of diffracted beam captured by ω scan

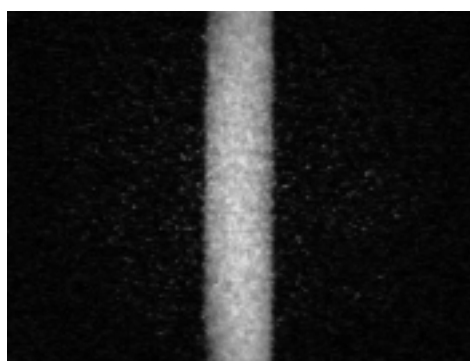

(b) X-ray diffraction image after adjusting orientation Si 440

Fig. 6.21 Axial adjustments

- (3) Setting the scanning range for the traverse stage
Scan the traverse stage while monitoring the image of the diffracted X-ray beam. The image of the diffracted beam disappears at the crystal edge. Stop scanning at this point. Do this at both ends of the crystal and register the scanning range.
- (4) Acquiring curve correction data when the sample is warped
If the sample is warped, the beam shifts from the diffraction condition with the scan of the traverse stage, and the X-ray image disappears. In such cases, acquire the step data for the curve correction and perform scanning with the correction. Here, the XRT-100 system supports only simple correction. If the sample is warped significantly at a curvature radius of 5 m or less, the correction cannot track the curvature and will generate an error.
- (5) Setting the variable slit
Set the precession, proceeding carefully to keep the slit from contacting the sample within the scanning range, and position as close to the sample as possible. Adjust the position of the right and left slits to control the aperture of the slit to avoid blocking the diffracted X-ray beam.
- (6) Attaching the film cassette
Insert an IP or an X-ray film into the cassette and attach the cassette to the cassette holder with the cassette bracket. Set the ψ angle along the axis of the diffraction angle by visual observation. Set the precession while taking steps to keep the cassette from contacting the third slit (S3) within the scanning range. Position as close as possible to the sample.
- (7) Developing exposed film
Set the conditions, including the number of scans, and expose the sample, then develop the film.
- (8) Devising imaging
Shift the ω angle so that it is slightly offset from the exact Bragg angle in imaging to increase the contrast of the topograph.
Vary the ω angle and take a series of topographs to acquire a topographical rocking curve containing angle information.
Scanning represents a type of averaging operation. Topography without scanning is also possible. Combined with the stepwise movement of the stage, this will produce a striped topograph.
Figs. 6.22 and 6.23 show examples of topographs.

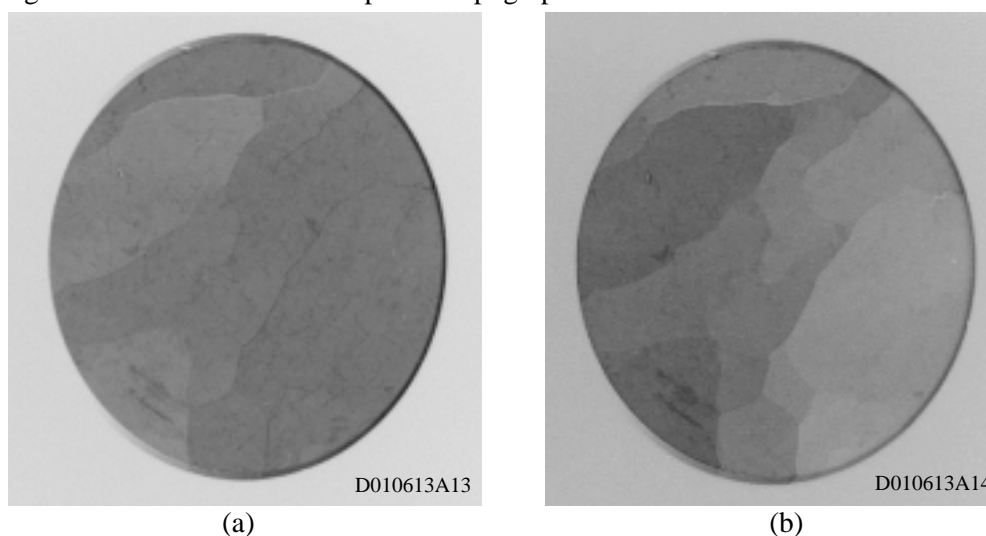


Fig. 6.22 Example of topography: effect of offset topography (negative)

Image (a) is taken at the exact Bragg angle, while Image (b) is taken slightly offset from the exact Bragg angle to the smaller angle. Offset imaging is more sensitive to changes in orientation. The diffraction vector is 440 .

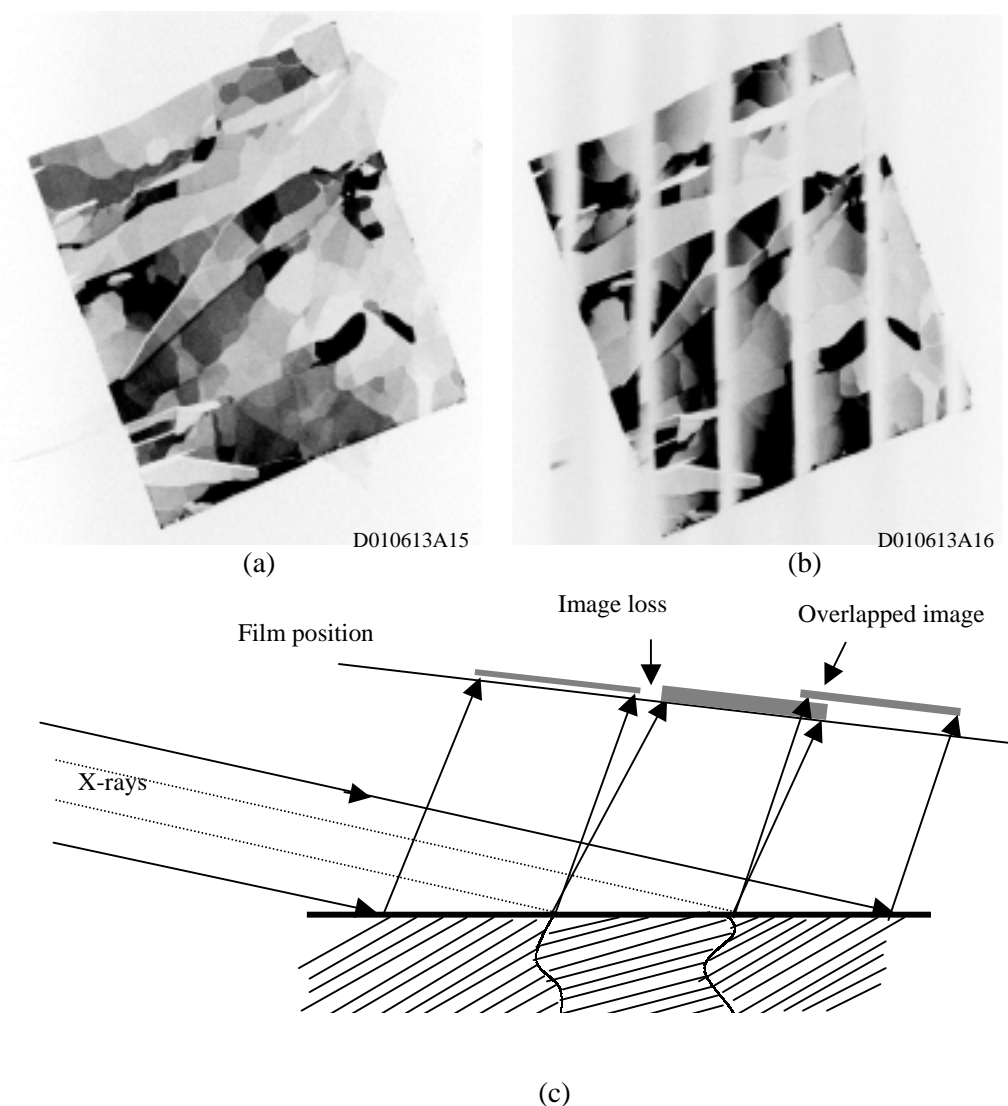


Fig. 6.23 Example of topography: striped topography (D010613A15,16) (negative)

Image (a) is a traverse topograph captured by the scanning technique. Image (b) is a topograph measured while the traverse stage is moved stepwise at constant intervals with repeated time exposures. The diffraction vector is 422 , and the topograph is captured with asymmetric reflection with an incident angle of approximately 14 degrees. The topographs clearly show the subgrain structure with orientation contrast. Image overlaps and image losses occur at the grain boundary due to changes along the axis of the diffracted beam as well as changes in the degree of diffraction conditions satisfied. Fig. 6.23 (c) schematically illustrates this behavior.

6.4.3 Scattered X-ray shielding plate for reflection topography

When measuring a small sample with a straight slit, attach the scattered X-ray shielding plate to prevent fogging caused by scattered X-rays. The plate extends the right piece of the third slit and shields against scattered X-rays.

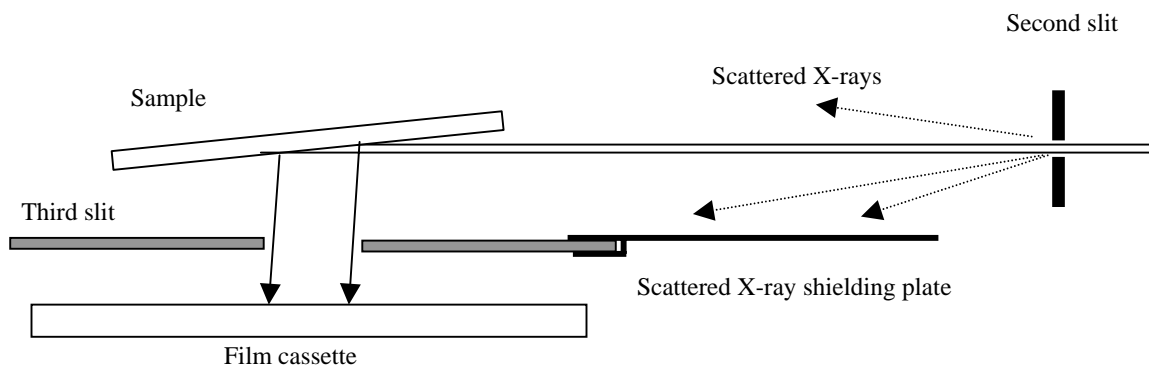


Fig. 6.24 Scattered X-ray shielding plate for reflection topography

6.5 Darkroom equipment

This section describes the darkroom equipment required for X-ray topography.

6.5.1 X-ray film and nuclear plates

(1) X-ray film

Fuji industrial X-ray film (manufactured by Fujifilm Corporation) may be used. Five types of films are available with Ix series numbers of #150, #100, #80, #50, and #25. Films with lower numbers use finer grains but have lower sensitivity. The #80 film is recommended as the standard film for X-ray topography. The size of the film is 10 inches by 12 inches (quartered). Twenty-five individual films are provided per box. Other sizes are also available: 3.1/3 inches × 12 inches and 14 inches × 17 inches (halved). Boxes with 50 films and 100 films are also available.

Kodak (US) and Agfa (Germany) manufacture equivalent products.

(2) Nuclear emulsion plates

Use nuclear plates manufactured by Ilford for topography. Ilford manufactures various types of emulsion, including L4 and G5. L4 incorporates the finest grains. The emulsion thicknesses best suited to topography are 25 μm and 50 μm . Plate dimensions are 1 inch × 3 inches. A standard box contains a dozen plates. Plates measuring 3 inches × 3 inches are also available.

(3) Using other films

FG film for electron microscope and medical film for mammography (high-resolution film for breast cancer examination) are single emulsion, high-resolution films, although with low sensitivity.

6.5.2 Developer

The specified X-ray film developer and fixer are listed below.

Developer: Rendol, 2 liters

Fixer: Renfix, 2 liters Both are manufactured by Fujifilm Corporation.

Dissolve in separate solutions of warm water to prepare developer and fixer solutions.

For the stopper solution, use a 1.5% aqueous solution of acetic acid (use Fuji acetic acid 11). Other solutions used include the Fuji QW (quick-washing) agent and Fuji Film Driwel Wetting agent.

Develop the nuclear emulsion plate with Kodak D-19 developer or ID-19 developer.

6.5.3 Darkroom equipment

Select the darkroom equipment from brochures provided by photo supply manufacturers (for example, Hansa by Hansa Co., Ltd. or King by Asanuma & Co., Ltd.).

Tools for development: Trays, tongs, jars for storing solutions, measuring beaker (2-liter), funnel, stirring rod, photo sponge, and thermometer

Darkroom facility: Ideally, the darkroom should contain a sink. A temporary space screened by blackout curtains or a space with basic darkroom facilities can be used as alternatives.

Darkroom equipment: Darkroom lamp (with various filters, including those for X-rays), darkroom timer, scissors, paper cutter (for cutting films),

- low-temperature thermostatic bath (for developing nuclear plates)
- Printing equipment: Enlarger, easel, focusing scope, monochromatic paper, developer and fixer solution for printing paper
- Scanner: The image can be transferred to a PC using a film scanner (e.g., Epson F-3200 film scanner for personal hobby use).
- Micrographs: It is also possible to take low-power micrographs. This is particularly useful for topographic observations with nuclear plates.

Also prepare tape for one-sided development (Scotch 845 Book tape or wide vinyl tape) of X-ray films.

7. Maintenance

7.1 Replacing the mylar films on the sample holder rings

- (1) Remove the broken mylar film.
Remove the adherend surface with the blade of a knife and clean the surface with a cleaner (e.g., benzene).
- (2) Prepare a cylinder with an inner diameter of 125 mm to 130 mm.
(Ideally, use an empty can (i). Open a hole to allow air to pass through.)
- (3) Stretch the mylar films (ii) tight at the bottom of the empty can and fasten with several dozen rubber bands (iii). Reinforce with Scotch tape.
- (4) Apply a thin coat of the adhesive provided with the system on the adherend mylar surface of the holder ring (iv), then place the mylar atop it. Place a weight (v) of approximately 5 kg to 6 kg on the mylar.
- (5) After an hour, the mylar film will have adhered completely. Remove excess mylar film with the blade of a knife.
- (6) After the mylar film is attached to each of the two holder rings, make approximately a dozen holes in the mylar film to allow air to pass through.

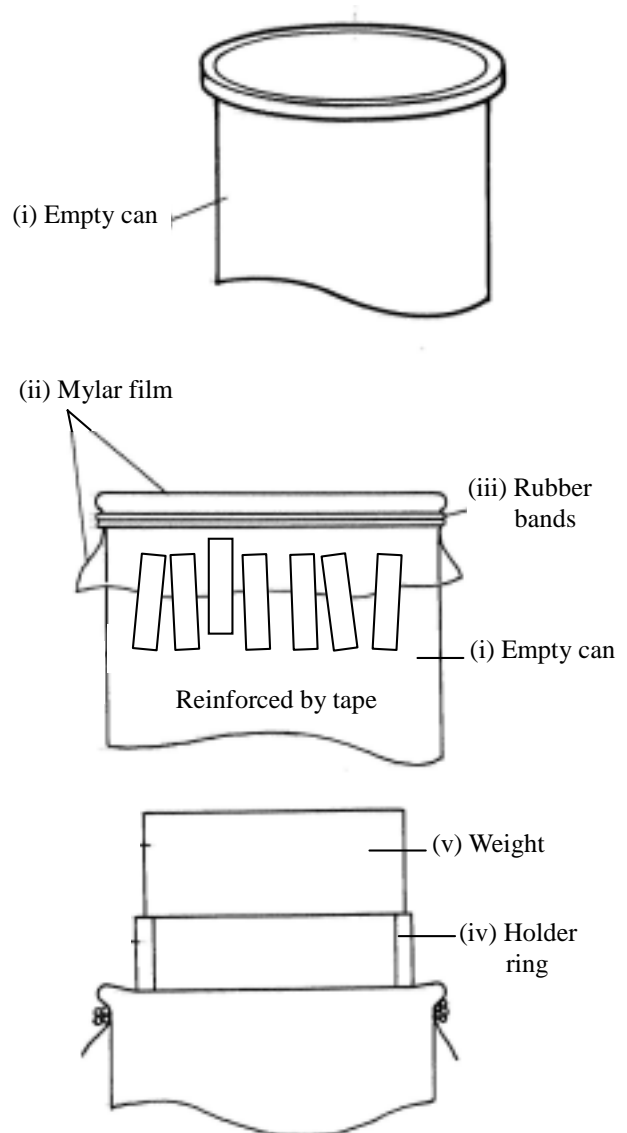


Fig. 7.1 Attaching mylar films

7.2 Installation

Position the components as indicated by the installation diagram. Place the mount for the main body on the floor. A bolt screw adjuster is provided to eliminate wobbling. Adjust the height of the beam center of the topogoniometer to 490 mm from the table. Use a level to ensure that the topogoniometer are level. Fine-adjust the height of the connecting section.

The radiation enclosure of the main body may be divided into halves. Reassemble after installing the mount for the main body.

Set the zero point as follows:

- (1) Set the 2θ rotary plate at the zero position;
- (2) Open the third slit wide to both sides;
- (3) Insert the 0.2-mm slit (for zero-setting) in the second slit holder and also in the fourth slit holder. Set the vertical position of the counter holder at zero;
- (4) Start the X-ray unit and align the camera with the two bottom screws, watching the fluorescent plate so that the beam that has passed the second slit runs through the fourth slit; and
- (5) Insert the 1-mm slit in the second slit holder and remove the height-limiting slit. Start the X-ray unit and observe the X-ray image with the fluorescent plate to check for an image cut in the vertical direction. If there is such a cut in the X-ray image, adjust the position of the X-ray shield cylinder until a uniform, full image is provided.

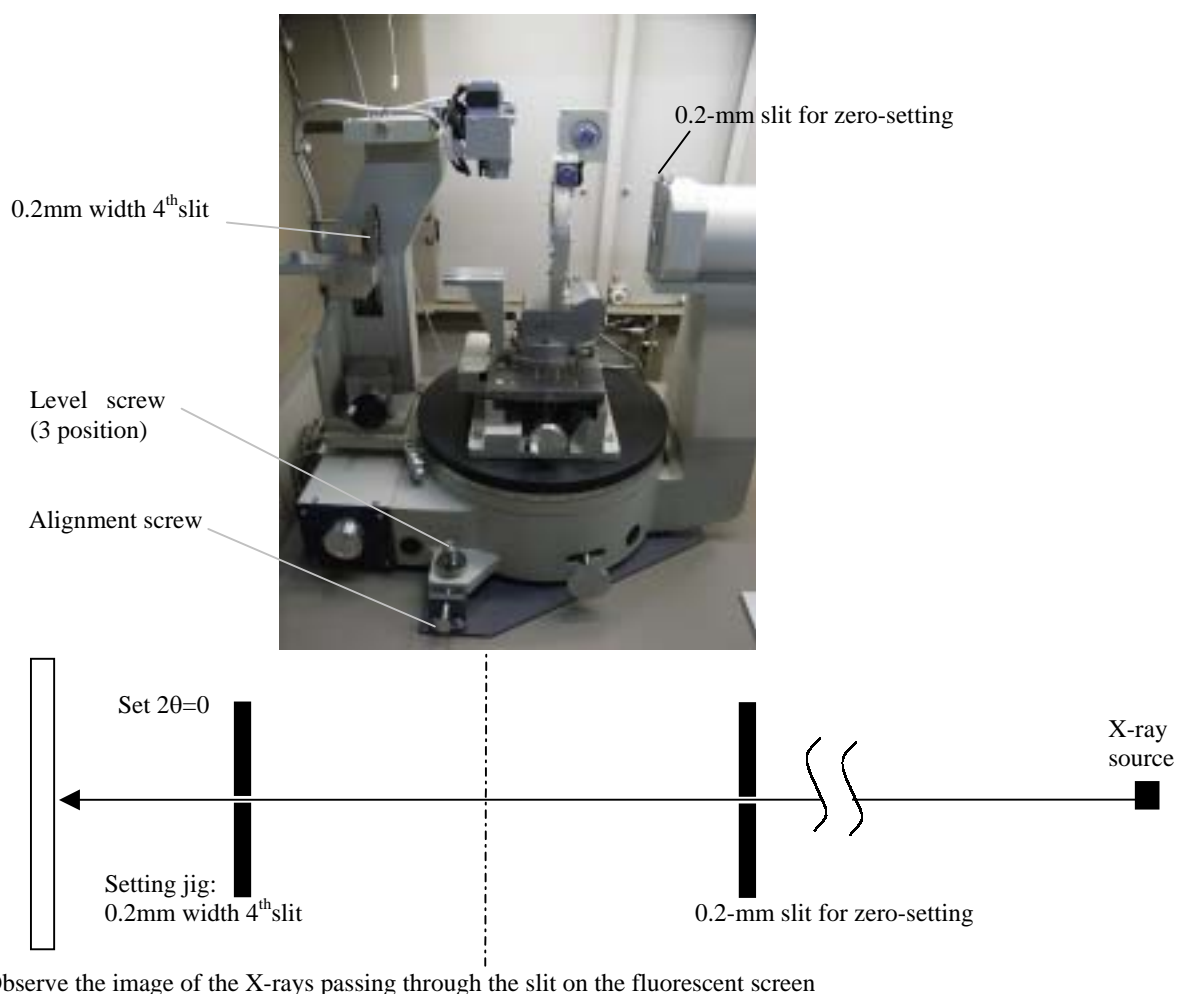


Fig. 7.2 Zero setting

Appendix

A Data

Table 1 2 B

(1) Si a = 0.5430nm

hkl	d (nm)	$2\theta_B (^\circ)$		
		AgK	MoK	CuK
111	0.3135	10.23	12.99	28.45
220	0.19198	16.75	21.29	47.31
311	0.1637	19.67	25.02	56.13
400	0.13575	23.78	30.29	69.13
331	0.12457	25.95	33.08	76.38
422	0.11084	29.23	37.31	88.04
511 333	0.10450	31.05	39.67	94.96
440	0.09599	33.88	43.36	$^\circ 106.72$
531	0.09178	35.48	45.45	114.11
$1/\mu$ (mm)		1.26	0.64	0.07

(2) Ge a = 0.5657nm

hkl	d (nm)	$2\theta_B (^\circ)$		
		AgK	MoK	CuK
111	0.32661	9.82	12.47	27.28
220	0.20000	16.08	20.43	45.30
311	0.17056	18.88	24.00	53.69
400	0.14143	22.81	29.04	66.00
331	0.12978	24.89	31.72	72.81
422	0.11547	28.03	35.77	83.68
511 333	0.10887	29.77	38.02	90.06
440	0.10000	32.48	41.54	100.75
531	0.0956	34.01	43.54	107.32
$1/\mu$ (mm)		0.055	0.029	0.025

(3) GaAs $a = 0.56534\text{nm}$

hkl	d (nm)	$2\theta_B (^\circ)$		
		AgK	MoK	CuK
111	0.32642	9.83	12.47	27.30
220	0.19989	16.09	20.44	45.33
311	0.17047	18.89	24.01	53.72
400	0.14135	22.82	29.06	66.04
331	0.12971	24.90	31.73	72.86
422	0.11541	28.05	35.79	83.74
511 333	0.10881	29.80	38.04	90.13
440	0.09995	32.50	41.57	100.83
531	0.09557	34.03	43.56	107.41
$1/\mu$ (mm)		0.051	0.027	0.023

Table 2 K Doublet

	Doublet		Natural breadth	
	d nm	d^2 nm ²	d nm	d^2 nm ²
AgK	0.000442	7.9×10^{-3}	0.000029	5.1×10^{-4}
MoK	0.000428	6.03×10^{-3}	0.000029	4.1×10^{-4}
CuK	0.000382	2.48×10^{-3}	0.000057	3.8×10^{-4}

B Guide for references related to topography

- A. Authier: “Dynamical Theory of X-ray Diffraction” revised edition, Oxford (2004) chapter 7
- D.K.Bowen & B.K.Tanner: High resolution X-ray Diffractometry and Topography, Taylor&Francis(1998)
- A. R. Lang: “International Tables for Crystallography Volume C” Part2, 2.7 Topography, KULWER ACADEMIC PUBLISHERS (1992)
- B.K.Tanner & D.K.Bowen ed.: “Characterization of Crystal Growth Defects by X- ray Method”, Plenum Press (1980)
- B.K.Tanner: “X-ray Diffraction Topography”, Pergamon Press (1976)