


Microscope Camera Quick Start


Micromanipulator Probe Station

January 27, 2021

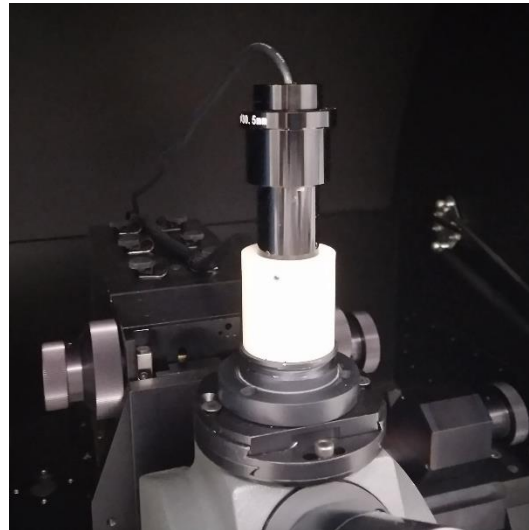


Camera Installation

The camera is installed at the top of the microscope; no adjustments should be necessary. Whatever you see looking through the eyepieces will be imaged by the camera.

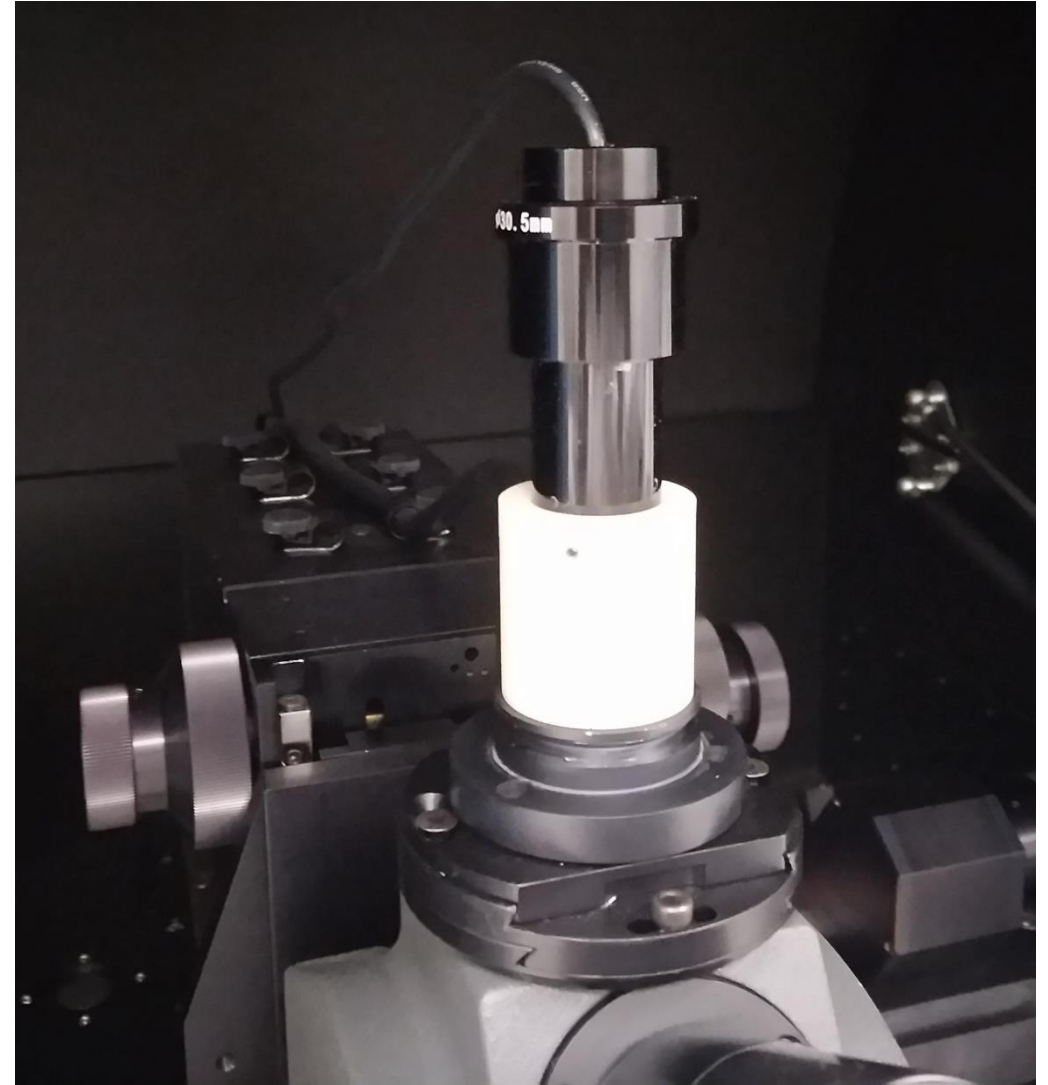

PC Desktop

Access the DinoCapture 2.0 Application by double clicking the icon

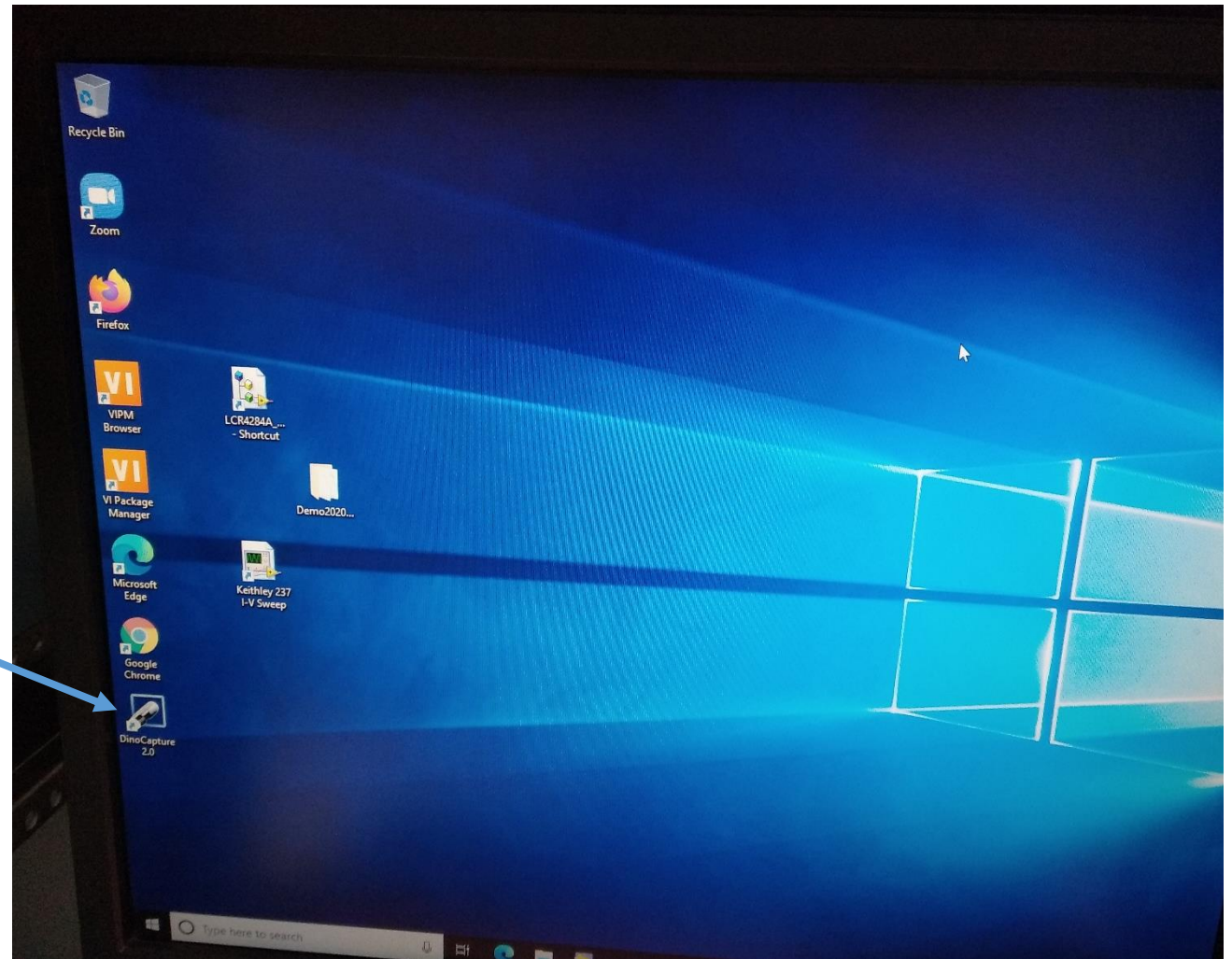


The camera interface is reasonably straightforward; you should be able to understand operation by experimentation...

Camera Controls

Button to take a picture


Live Image

Camera properties dialog

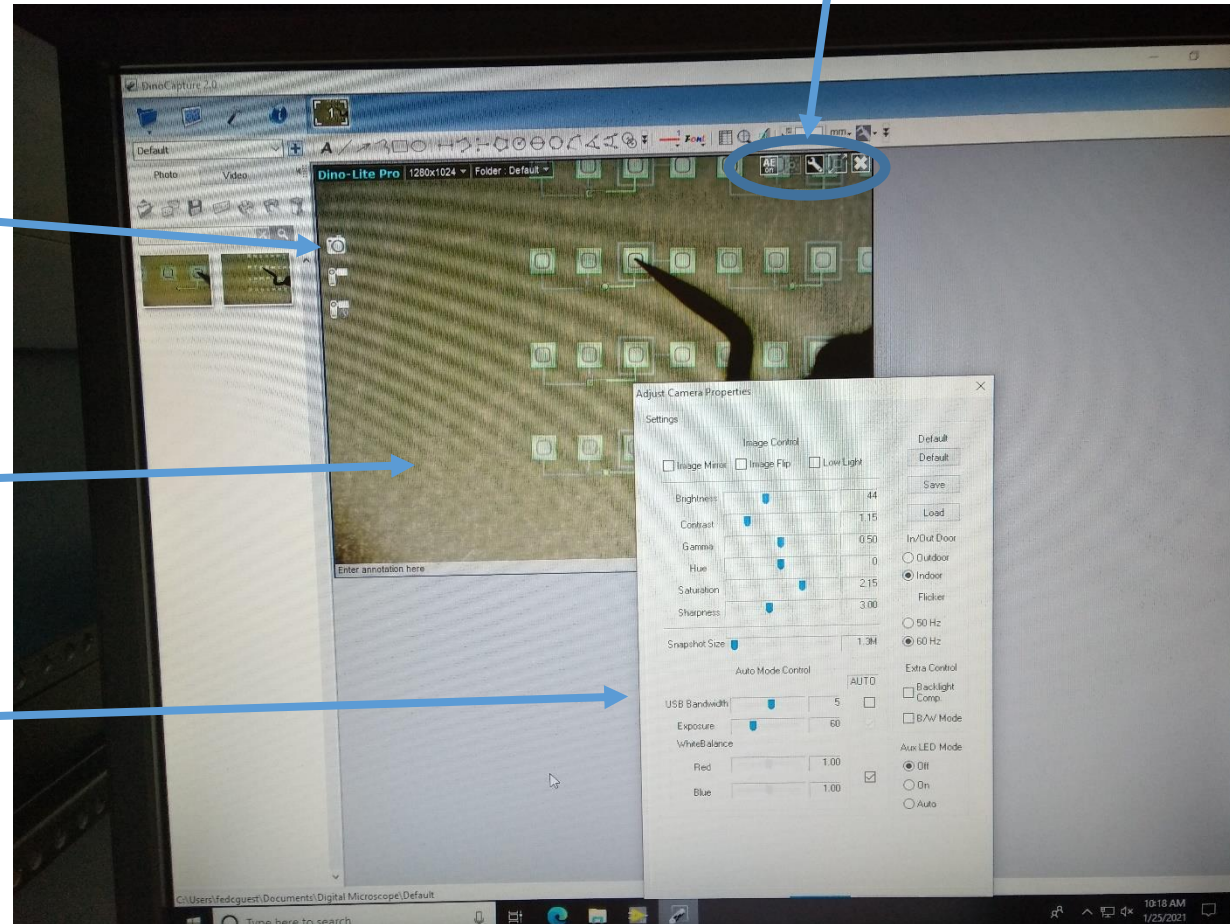

You can experiment with the camera properties to see how each one affects the image. Usually the defaults work well, however.

If image shows bands of noise, adjusting the USB bandwidth to a lower number may be helpful. The camera response will slow significantly however.

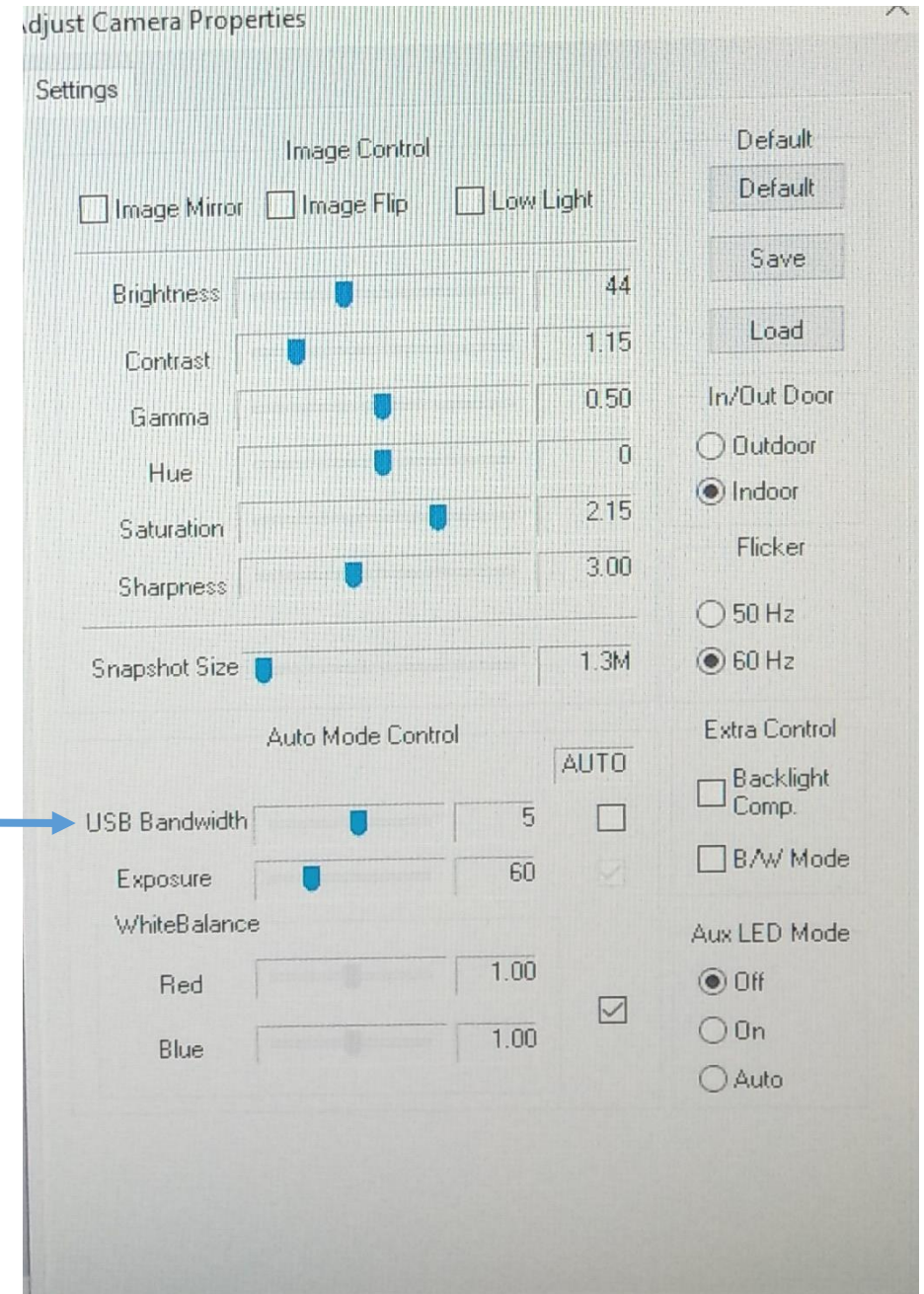

Sample Image Captures

